



10. Wall mount
11. Magnet
12. Table stand
13. Battery door

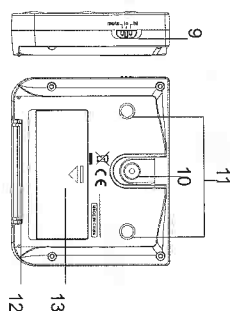

## LCD DISPLAY

1. Timer mode indicator
2. Timer1 window
3. Timer2 window
4. Hour
5. Minute
6. Second
7. Count down indicator
8. Count up indicator

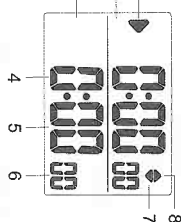


## COUNT DOWN TIME MEMORY SETTING

1. In Timer1/Timer2 mode, press and hold "M" key for 2 seconds to enter count down time memory setting.
2. Press "HR", "MIN", "SEC" key to set the Hour, Minute and Second separately.
3. Press "M" key again to confirm and exit the setting, or let it exit automatically 30 seconds later without pressing any key.

### Note:

The clock will store the time value as memory time.

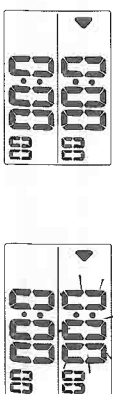

Count Time Mode Count Down Time Memory Setting

## USING COUNT DOWN TIME MEMORY FUNCTION

1. Set the count down time memory as described in the previous section.
2. Press "S/S" key once to start the countdown function.
3. Press "S/S" key again to stop the countdown function.
4. The alarm will ring when the time count down to 0, press "S/S" key to stop the alarm and return to the pre-set memory time.

### Note:

1. During counting time mode, the green light indicator will keep flashing.

2. The clock will turn to count up time mode automatically when the countdown time to 0. With the alarm ringing synchronized.
3. The alarm sound will last for 1 minute, with red light indicator flashes.

## USING COUNT DOWN TIME FUNCTION


1. In Timer1/Timer2 mode, set count down time by pressing "HR", "MIN", "SEC" key to set the Hour, minute and second separately.
2. Press "S/S" key to start the countdown function.
3. Press "S/S" key again to stop the countdown function.
4. The alarm will ring when the time count down to 0, press "S/S" key to stop the alarm and return to the pre-set count down time.

## USING COUNT UP TIME FUNCTION

1. When Timer1/Timer2 displays 0:00 00, press "S/S" key once to start count up time function.
2. Press "S/S" key again to stop the count up function.

### Note:

The timer will count up to 99 hour 59 minute 59 second unless press "S/S" key to stop it.

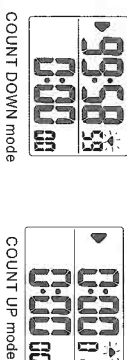

COUNT DOWN mode

COUNT UP mode

## HOW TO CHANGE THE COUNT DOWN TIME MEMORY

1. Press and hold "M" key for 2 seconds until the digits flashes, and then press "HR", "MIN", "SEC" key to set the desired value.
2. Press "M" key again to confirm and exit the setting.


## HOW TO CLEAR THE DATA

Set the count down time and count down time memory as described in the previous sections, press "HR" and "MIN" key at the same time to clear all the data to 0 after stop time counting.

## HOW TO PLACE THE CLOCK

1. This clock has both desktop and wall-mounting structures.

Place the screw on the desired wall and hang the clock by the wall mount in the back of it or just simply place it on the desktop by its table stand. Furthermore, you can use the magnet in the back of the clock to attach to the metal surface.

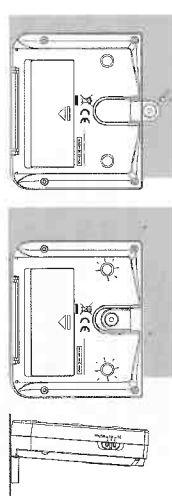

## BATTERY REPLACEMENT

When the LCD becomes dim, replace with two new AAA size batteries at once. Pay attention to the battery polarity.

### Note:

Attention! Please dispose of used unit or batteries in an ecologically safe manner.

