

Histamine ELISA Test Kits

The AgraQuant® Histamine and AgraQuant® Histamine Rapid test kits are competitive enzyme-linked immunosorbent assays (ELISA) for the sensitive and quantitative analysis of histamine in a variety of food products.



AgraQuant® Histamine

The AgraQuant® Histamine ELISA Test kit combines an acylation process with a competitive ELISA assay designed for the quantification of histamine in fish meal, fresh fish, milk, cheese, sausage and wine.

By use of the acylation reagent, any histamine present in the sample is derivatized into N-acylhistamine, which can subsequently be quantified with the ELISA method.

The sample preparation and extraction procedure have been optimized for the determination of histamine in some of the most common matrices.

Matrix	Quantitation Range	
Fish meal	100 - 10000 ppm	
Fish, cheese, sausage	2.5 – 250 ppm	
Milk	0.1 - 100 ppm	
Wine, champagne	0.25 - 25 ppm	

AgraQuant® Histamine Rapid

The AgraQuant® Histamine Rapid ELISA Test kit employs a competitive ELISA assay for the quantitative determination of histamine in fresh, canned and salted fish, as well as in fish meal and fish in oil.

The workflow of this assay has been optimized for the analysis of histamine in fish and therefore allows for a much faster assay procedure.

With only 15 minutes necessary for the sample preparation and a total incubation time of 20 min for the ELISA assay, the results can be obtained as fast as in 35 minutes.

Matrix	Quantitation Range
Fish meal	3 – 300 ppm
Fresh and other fish	3 – 300 ppm

PRODUCT FEATURES

Highly specific and sensitive

 Complies with regulatory limits of histamine in food

Easy

Simple sample preparation and ready to use reagents

Cost-effective

Minimizes waste and maximizes value



AgraQuant® Histamine Test Kits

Please read the instructions on the package insert completely before performing any test.

3-Step Procedure



Extraction

Extract your sample with distilled water by following the simple step-by-step instructions given in the package insert. There are clear descriptions for each matrix.



Acylation

By use of the acylation reagent, histamine in the food extracts is quantitatively derivatized into N-acyl-histamine. All required materials are supplied within the kit.



ELISA Assay

The competitive Histamine ELISA kit uses the microtiter plate format. Histamine is coated to the solid phase of the microtiter plate. Acylated histamine and solid-phase-coated histamine compete for the antiserum binding sites. When the system is in equilibrium, free antigen and free antigen-antiserum complexes are removed by the washing step. The antibody bound to the solid-phase coated histamine is then detected by anti-goat IgG conjugated with peroxidase (enzyme conjugate). An enzyme substrate is added and blue color develops. A stop solution is then added to change the color from blue to yellow. The microwells are measured optically using a microwell reader with a reading filter of 450 nm and a differential filter between 620 nm to 650 nm.

Ordering Information - Romer Labs Histamine Test Kits

Product	Format	Item Number	Quantitation Range
AgraQuant® Histamine	96 wells	10002017	see front page*
AgraQuant® Histamine Rapid	48 wells	10002018	see front page*

^{*} matrix-dependent

Romer Labs Inc. - America · Newark, DE, USA · T +1 302 781 6400 · F +1 302 781 6378 · E office@romerlabs.com
Romer Labs Diagnostic GmbH - Europe · Tulln, Austria · T +43 2272 615 33 · F +43 2272 615 33 13177 · E office-europe@romerlabs.com
Romer Labs Singapore Pte. Ltd. - Asia · Singapore · T +65 6631 8018 · F +65 6275 5584 · E salesasia@romerlabs.com