Compact Dry EC medium for E.coli and coliforms/Medium für E.coli und Coliforme/milieu pour E.coli et coliformes/medio para E.coli y coliformes/ medio per E.coli e coliformi / meio para E.coli e coliformes

40 plates/Platten/plaques/placas/lastre/placas

ID-No. 1 000 168

240 plates/Platten/plaques/placas/lastre/placas

ID-No. 1 000 169

920 plates/Platten/plaques/placas/lastre/placas

ID-No. 1 002 878

English

Deutsch

Français

Compact Dry EC is a ready to use, chromogenic plate for the enumeration of E.coli and coliforms

Sample pretreatment

Viable count in water or liquid foodstuff

Drop 1 ml of specimen (dilute if necessary) on the middle of the Compact Dry plate.

Viable count in solid foodstuff

Add buffer solution to the sample and homogenize by stomacher®. Drop 1 ml of specimen (dilute if necessary) on the middle of the dry sheet of the Compact Dry plate.

Viable count in swab test specimen

Use the swab to wipe the surface, put into the device with wiping solution. Drop 1 ml of wiping solution (dilute if necessary) on the middle of the Compact Dry plate. It is recommended to use "Swab for Compact Dry" offered by HyServe Id-No. 1 002 952/3 (40/240 pieces)

Test instructions

1. Open the cap and drop 1 ml of specimen on the middle of the Compact Dry plate.
2. Specimen diffuses automatically and evenly into the sheet and transforms the dried sheet into a gel within seconds.
3. Put the cap again on the plate and write the information needed on the memorandum section.
4. Turn over the capped plate and put in the incubator.
5. After incubation count the number of colored colonies underneath the plate. White paper placed under the plate helps to count the colonies.

Incubation time 24 ± 2 hours

Incubation temperature 37 ± 1°C

Interpretation of the results

E.coli forms blue colonies, red and blue colonies together are the total coliform group count.

Storage and shelf life

Keep at room temperature (+ 1 - 30 °C). Total shelf life 24 months after manufacturing.

Notes

- E. coli 0157 forms pink/red purple colonies.
- High concentrations on plates will cause the entire growth area to become white/pink. In this case dilute the sample.
- After use please follow the current disposal regulations.
- The growth area is 20 cm². The back of the plate has a grid carved of 1 cm x 1 cm to make the colony counting easier. In case of any difficulties to count colonies due to large number of colonies grown, the bacteria count can be obtained by multiplying 20 by an average number of colonies per grid counted from several grids. However, dilutions are recommended.
- Compact Dry plates are produced at an ISO 9001 certified site.
- AOAC approved; certificate No. 110402 MicroVal approval No.0805-004LR according to ISO 4832 (2006) and No.0805-005LR according to ISO 16649-2-2:2001
- ISO EN 16140:2003
- NordVal approval No. 036 and 037

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