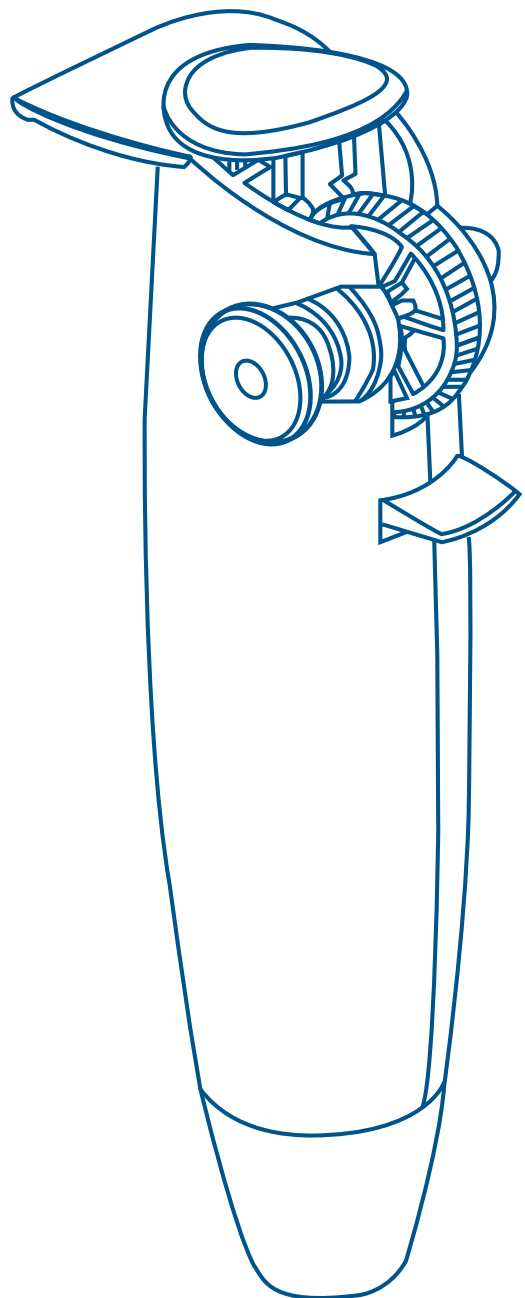
Instruction manual

VWR® Smoothie

North American Catalogue Number: 75856.440

European Catalogue Number: 612-4546

CE

VWR [®]
COLLECTION

Table of contents

EN Instruction Manual 1 - 6	DE Bedienungsanleitung 7 - 12	IT Manuale d'istruzioni 13 - 18
FR Manuel d'instructions 19 - 24	ES Manual de instrucciones 25 - 30	

United States

Manufactured for:

VWR International, LLC
100 Matsonford rd
Radnor, PA 19087
800 932 5000
us.vwr.com

Europe

Address of legal manufacturer:

VWR International bvba
Researchpark Haasrode 2020
Geldenaaksebaan 464
B-3001 Leuven
+ 32 16 385011
be.vwr.com

Contents

OPERATOR INSTRUCTIONS

4

INTRODUCTION

4

FITTING THE PIPETTE

4

**CHOOSING THE CORRECT THUMB
WHEEL SETTING**

4

FILLING THE PIPETTE

5

DISPENSING

5

'BLOW-OUT' PIPETTES

5

Operator Instructions

Introduction

The Smoothie pipette filler may be used with glass or plastic pipettes to facilitate the pipetting of a wide range of liquids. Pipettes within the volume range of 1 ml to 25 ml can be used. When correctly used, the liquid should only come into contact with the pipette.

Operating Instructions

1. Fitting the pipette

Carefully attach the pipette by inserting it into the silicone holder inside the nose cone. Do not use force.

2. Choosing the correct thumb wheel setting

The Smoothie is equipped with a rotating thumb wheel which advances the piston up and down the barrel for aspiration and dispensing. The unique thumb wheel can be set at two different dispense resolutions for maximum accuracy depending on the volume of liquid being dispensed. Fast filling for volumes of 10 ml to 25 ml is possible when the thumb wheel is positioned to the right. This is the default setting. Whenever you are working with pipettes smaller than 10 ml we recommend that you choose the fine setting – whereby the thumb wheel is positioned to the left.

To change the thumb wheel setting the piston must be in its lowest position (image 1A). To change from the default setting to the fine setting use the thumb of your right hand to push the thumb wheel by pressing the thumb wheel button all the way in to the left (image 1B). With your left hand rotate the fixing button until the gear locator is locked in its respective groove (image 1C), thus holding the thumb wheel in its new position. The thumb wheel should now be positioned to the left. The pipette is now ready for work in its fine setting as recommended for volumes less than 10 ml.

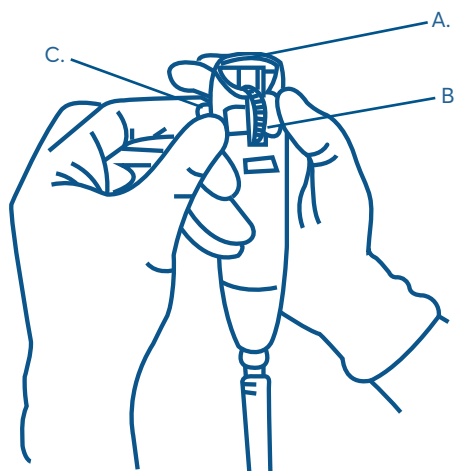
Image 1.

Choose the correct thumb wheel

- A. Piston in low position
- B. Push
- C. Rotate the fixing button

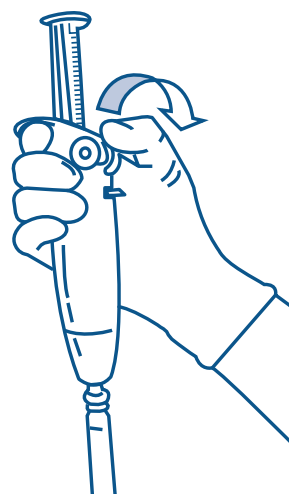
Image 2.

Rotate the thumb wheel to fill the pipette.

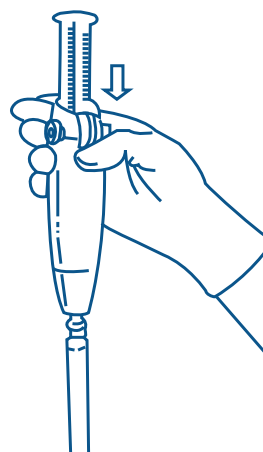
Image 3.

Press the dispense lever down to empty the pipette by gravity.

Operator Instructions

3. Filling the pipette

Place the tip of the pipette into the liquid to be aspirated. With your thumb rotate the thumb wheel (as shown in image 2) allowing the liquid to fill the pipette so that the meniscus is level with the required mark.

Note: Do not allow liquid to enter the Smoothie.

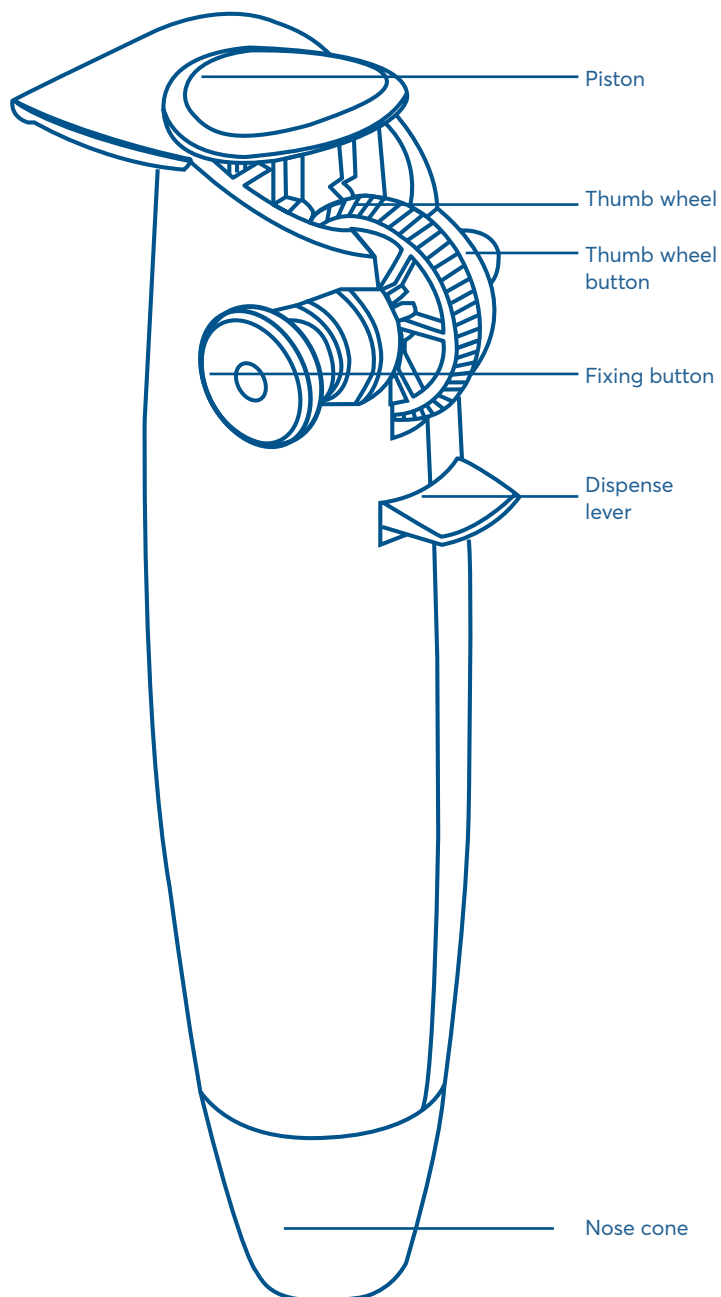
4. Dispensing

Hold the receiving vessel in an inclined position, placing the pipette tip against the receiving vessel's wall. Press the dispense lever down to empty the pipette by gravity (as shown in image 3).

5. 'Blow-out' Pipettes

As soon as the meniscus in the pipette comes to a standstill, rotate the thumb wheel upwards to expel the last few drops. To ensure the complete volume is dispensed, wipe the pipette tip upwards a few millimetres along the inside wall of the receiving vessel.

Using your free hand push the piston to its lowest position. The Smoothie is ready for a new pipetting operation.

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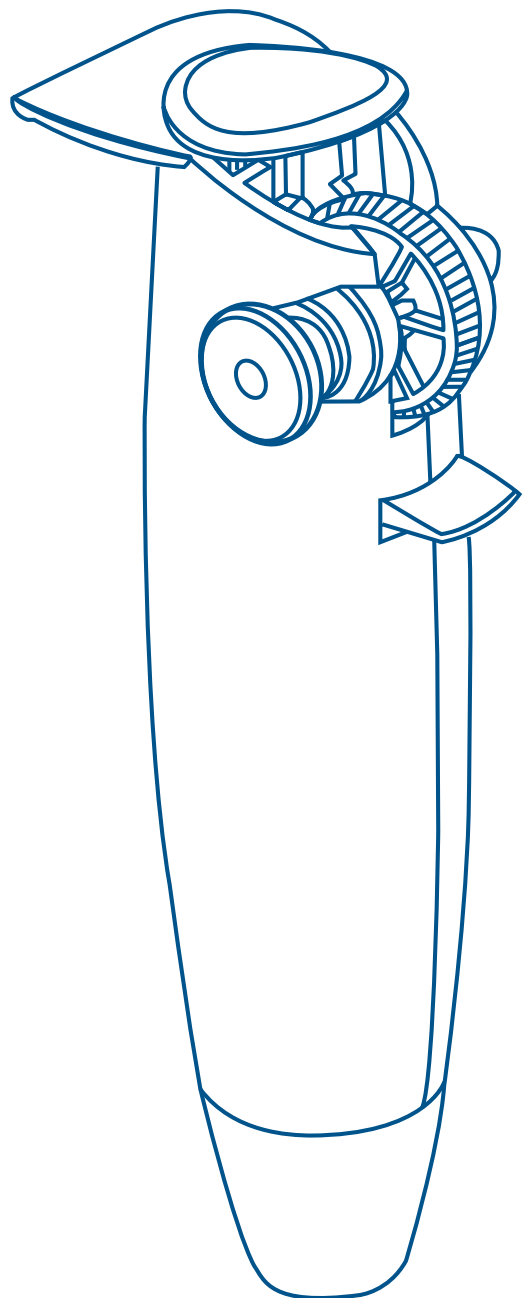
IRELAND / NORTHERN IRELAND

VWR International Ltd / VWR
International (Northern Ireland) Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: +353 (0) 1 88 22 222
sales.ie@vwr.com

ITALY

VWR International S.r.l.
Via San Giusto 85
20153 Milano (MI)
Tel.: +39 02 3320311
info.it@vwr.com

THE NETHERLANDS

VWR International B.V.
Postbus 8198
1005 AD Amsterdam
Tel.: +31 (0) 20 4808 400
info.nl@vwr.com

NORWAY

VWR International AS
Brynsalleen 4,
0667 Oslo
Tel.: +47 22 90 00 00
info.no@vwr.com

POLAND

VWR International Sp. z o.o.
Limbowa 5
80-175 Gdansk
Tel.: +48 58 32 38 200
info.pl@vwr.com

PORTUGAL

VWR International - Material de
Laboratório, Lda
Centro Empresarial de Alfragide
Rua da Indústria, nº 6
2610-088 Amadora
Tel.: +351 21 3600 770
info.pt@vwr.com

SPAIN

VWR International Eurolab S.L.
C/ Tecnología 5-17
A-7 Llinars Park
08450 - Llinars del Vallès
Barcelona
Tel.: +34 902 222 897
info.es@vwr.com

SWEDEN

VWR International AB
Fagerstagatan 18a
163 94 Stockholm
Tel.: +46 (0) 8 621 34 00
kundservice.se@vwr.com

SWITZERLAND

VWR International GmbH
Lerzenstrasse 16/18
8953 Dietikon
Tel.: +41 (0) 44 745 13 13
info.ch@vwr.com

UK

VWR International Ltd
Customer Service Centre
Hunter Boulevard - Magna Park
Lutterworth
Leicestershire
LE17 4XN
Tel.: +44 (0) 800 22 33 44
uksales@vwr.com

CHINA

VWR (Shanghai) Co., Ltd
Bld.No.1, No.3728 Jinke Rd,
Pudong New District
Shanghai, 201203- China
Tel.: 400 821 8006
Email: info_china@vwr.com

INDIA

VWR Lab Products Private Limited
No.139, BDA Industrial Suburb,
6th Main, Tumkur Road, Peenya Post,
Bangalore, India - 560058
Tel.: +91 80 28078400
vwr_india@vwr.com

MIDDLE EAST & AFRICA

VWR International FZ-LLC
DSP Laboratory Complex
125, Floor 01
Dubai, United Arab Emirates
Tel.: +971 4 5573271
info.mea@vwr.com

SINGAPORE

VWR Singapore Pte Ltd
18 Gul Drive
Singapore 629468
Tel.: +65 6505 0760
sales.sg@vwr.com

GO TO VWR.COM FOR THE LATEST NEWS, SPECIAL OFFERS AND DETAILS FROM YOUR LOCAL VWR SUPPORT TEAM