


Application note for the flexible Universal Drain Tube for OPTIFIX® SAFETY, SAFETY S and OPTIMAT® 3 Art.-No. 101.104

Instead of the discharge tube set, the universal drain tube is screwed onto the thread at the valve block.

In order to deaerate the device and the universal drain tube, the finger protection cap which encloses the dosing piston is raised and pressed down a number of times so that no air bubbles are visible in the dosing cylinder or in the universal drain tube.

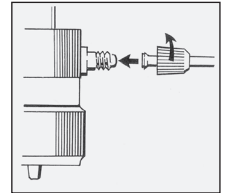

Please note that even when using the universal drain tube you have to set the desired dosing volume at the volume setting system.

Insert the universal drain tube into the dosing vessel, pull the finger protection cap up and dose the desired amount of liquid by pressing the finger protection cap down.



All data are subject to change • Copyright Poulten & Graf GmbH, Karl-Carstens-Str. 10, D-97877 Wertheim - 2017

Distribué par :

Z.A de Gesvrine – 4 rue Képler – B.P.4125
44241 La Chapelle-sur-Erdre Cedex – France
t. : +33 (0)2 40 93 53 53 | f. : +33 (0)2 40 93 41 00
commercial@humeau.com

 Laboratoires
Humeau

w w w . h u m e a u . c o m