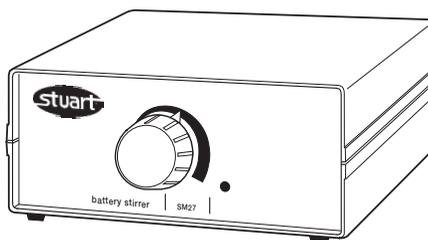
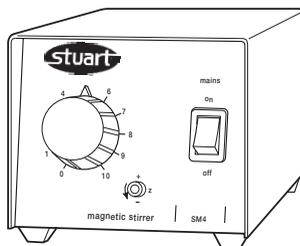
Magnetic Stirrer SM4

Portable Magnetic Stirrer SM27

Instructions for use
Instructions d'utilisation
Istruzioni per uso
Instrucciones de utilización
Bedienungsanleitung

English

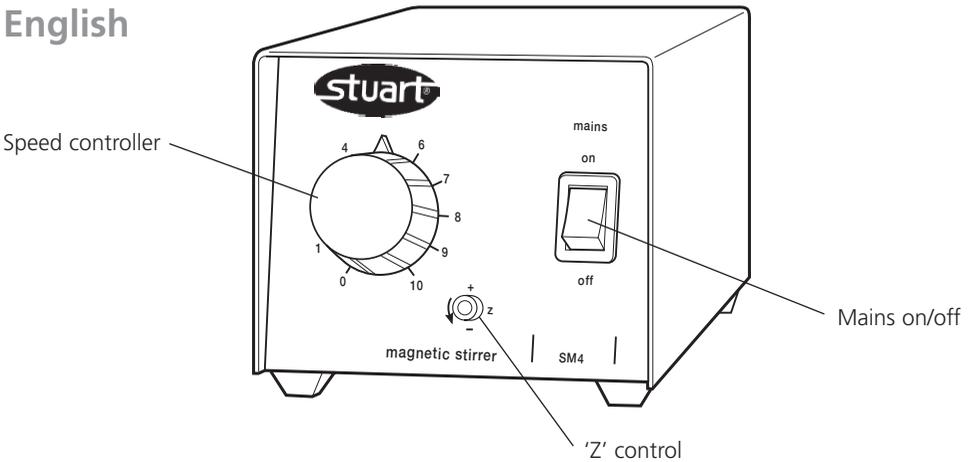
Figure 1: SM4 magnetic stirrer

SM4 MAGNETIC STIRRER

Thank you for purchasing this Stuart product. To get the best performance from the equipment, and for your own safety, please read these instructions carefully before use.

This equipment is designed to operate under the following conditions: -

- ❖ For indoor use only
- ❖ Use in a well ventilated area
- ❖ Ambient temperature +5°C to +40°C
- ❖ Altitude to 2000m
- ❖ Relative humidity not exceeding 80%
- ❖ Mains supply fluctuation not >10%
- ❖ Overvoltage category II IEC60364-4-443
- ❖ Pollution degree 2 IEC664
- ❖ Use with a minimum distance all around of 200mm from walls or other items

Electrical Installation

THIS EQUIPMENT MUST BE EARTHED

Before connection please ensure that the line supply corresponds to that stated on the rating label. The SM4 model requires a supply rated at 220-240Vac, 50Hz single phase.

There is an IEC socket at the rear on the left hand side of the instrument for connection to the mains supply, (see figure 1). The unit is supplied with two mains leads fitted with IEC plugs for connection to the instrument. One has a U.K. 3 pin plug and the other has a 2 pin "Shuko" plug for connection to the mains supply. Choose the lead appropriate for your electrical installation and discard the other.

Should neither lead be suitable you should obtain a moulded lead locally. If this is not possible, take the lead with the U.K. plug and replace the plug with a suitable alternative.

Should the mains lead require replacement a cable of 1mm² of harmonised code H05V V-F connected to an IEC 320 plug should be used.

N.B. The UK mains lead is protected by a 10A fuse mounted in the plug top.

The mains lead should be connected to the instrument BEFORE connection to the mains supply.

IF IN DOUBT CONSULT A QUALIFIED ELECTRICIAN

Safety Advice Before Use

- The unit should be carried using both hands with fingers under each side frame.
- Never move or carry the unit with containers on the top plate or while the unit is connected to the mains electricity supply.
- There is a danger of liquid spillage and splashing if containers are over filled and stirred at high speed. Always build the stirrer speed up slowly and never stir more rapidly than is necessary.

SM4 General Description

The SM4 is a mini magnetic stirrer ideal for general laboratory use and is shown in figure 1.

Operation

Position the apparatus on a flat surface.

Ensure the speed control is at the 'OFF' position. Place a magnetic follower in the vessel and position it centrally on the plate. Switch the unit 'ON' and turn the speed controller slowly to the right until the follower is rotating to the desired speed.

The lowest speed setting can be adjusted using the 'Z' Control. Turn the speed controller knob to number 1 and adjust the 'Z' Control until the follower is rotating at the desired speed. Every time the speed controller is set to 1, this speed will be observed.

Maintenance & Servicing

 **WARNING: Ensure the unit is disconnected from the mains electricity supply before attempting any cleaning or servicing.**

Cleaning

Periodically clean the instrument using a damp cloth and mild detergent solution. Do not use harsh or abrasive cleaning agents.

Servicing

This unit does NOT require routine servicing.

Repairs

Only spare parts supplied or specified by Bibby Sterilin Ltd. or its agents should be used. Fitting of non-approved parts may affect the performance and safety features designed into the instrument.

Any repairs or replacement of parts MUST be undertaken by suitably qualified personnel.

If in any doubt, please contact the Technical Dept. of Bibby Sterilin or the point of sale.

Bibby Sterilin Ltd.

Stone

Staffordshire,

ST15 0SA

United Kingdom

Tel: +44 (0) 1785 812121

Fax: +44 (0) 1785 813748

Technical Specification

Speed, range	80 to 1300rpm
Overall dimensions (w x d x h)	90 x 160 x 90mm
Net weight	1kg
Electrical supply	230V, 50Hz, 50W

Warranty

Bibby Sterilin Ltd. warrants this instrument to be free from defects in material and workmanship, when used under normal laboratory conditions, for a period of two (2) years. In the event of a justified claim Bibby Sterilin will replace any defective component or replace the unit free of charge.

This warranty does NOT apply if damage is caused by fire, accident, misuse, neglect, incorrect adjustment or repair, damage caused by incorrect installation, adaptation, modification, fitting of non-approved parts or repair by unauthorised personnel.

SM27 PORTABLE MAGNETIC STIRRER

Thank you for purchasing this Stuart product. To get the best performance from the equipment, and for your own safety, please read these instructions carefully before use.

This equipment is designed to operate under the following conditions: -

- ❖ For indoor use only
- ❖ Use in a well ventilated area
- ❖ Ambient temperature +5°C to +40°C
- ❖ Altitude to 2000m
- ❖ Relative humidity not exceeding 80%
- ❖ Mains supply fluctuation not >10%
- ❖ Over voltage category II IEC60364-4-443
- ❖ Pollution degree 2 IEC664
- ❖ Use with a minimum distance all around of 200mm from walls or other items

If the equipment is not used in the manner described in this manual the protection provided by the equipment may be impaired.

Battery information

FOUR batteries, 1.5V, size D, IEC no. LR20

The batteries are fitted as shown in figure 2, with the positive terminals facing the rear of the instrument:

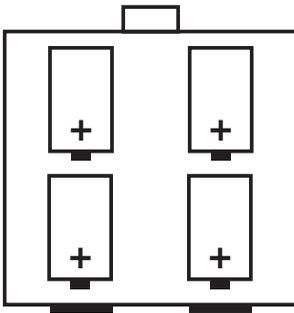
Figure 2: Battery orientation

Standard Alkaline Magnesium cells are recommended, but Zinc Chloride or rechargeable Nickel Cadmium batteries can also be used. Rechargeable batteries MUST be removed from the stirrer before recharging.

Main adapters

The SM27 stirrer can also be operated from an external 230Vac power supply PROVIDED THAT THE APPROPRIATE MAINS ADAPTER IS USED. Suitable adapters are detailed under 'Spares & Accessories'.

Adapters can be used with or without the batteries in place.

WARNING: The SM27 MUST NOT be connected directly to the mains electricity supply.

IF IN DOUBT CONSULT A QUALIFIED ELECTRICIAN

Safety Advice Before Use

- Never move or carry the unit with containers on the top plate.
- Do not leave discharged batteries in the unit as these may leak and cause corrosion.
- Remove rechargeable batteries before recharging
- Use only an appropriate adapter if utilising a power source other than batteries.

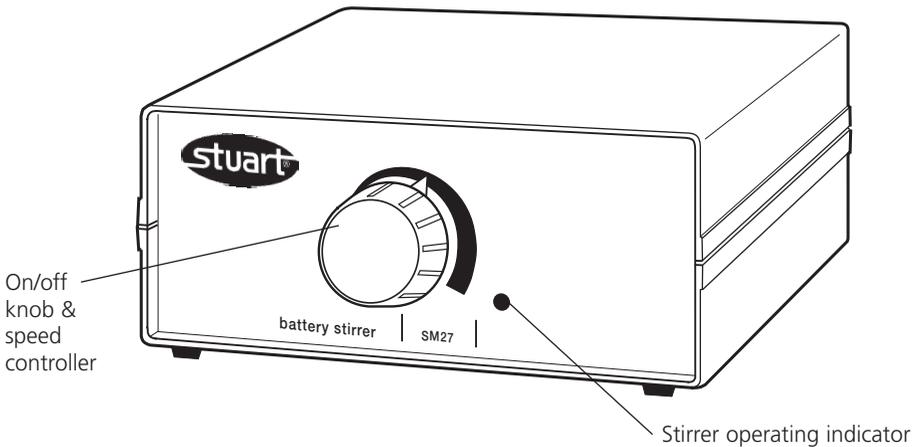
Figure 3: SM27 front view

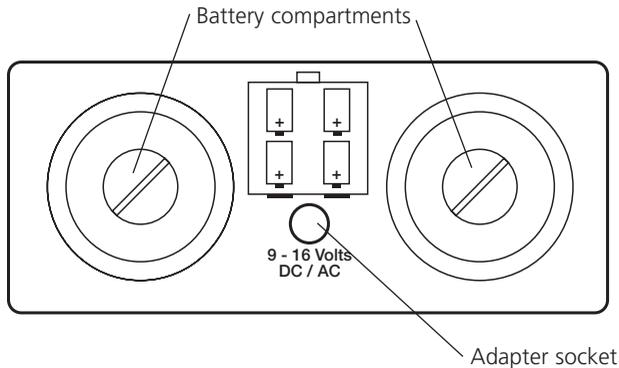
Figure 4: SM27 rear view

SM27 General Description

The SM27 stirrer shown in figure 3 can be battery powered or used on mains supply. It is ideal for use inside incubators, glove boxes or in the field.

Operation

Stirring

Place the solution to be stirred together with a PTFE follower in a suitable vessel and place in the centre of the unit. Switch the stirrer 'ON' and adjust the speed of stirring using the speed control knob (figure 3). When the unit is operating, the LED will be lit (figure 3).

Replacing the batteries

When the batteries are running low, the stirrer operating indicator (figure 3) will dim. To replace the batteries, open the battery compartments (figure 4) using a coin. Remove used batteries and dispose of safely. Replace new batteries with positive terminals facing the rear as shown in figures 2 & 4.

Using an adapter

Optional external power supply adapters are available for 230Vac and a car cigarette socket (see accessories overleaf).

Connect the stirrer to the adapter using the low voltage socket at the back of the unit (figure 4) before connecting the power supply.

Maintenance and Servicing

WARNING: Ensure the batteries are removed and /or the unit is disconnected from the mains adapter before attempting maintenance or servicing.

Cleaning

Periodically clean the instrument using a damp cloth and mild detergent solution. Do not use harsh abrasives or abrasive cleaning agents. Avoid strong acids and solvents on the ABS case.

Spares and Accessories

The following spares and accessories are available from Bibby Sterilin or your laboratory dealer:

Description	Catalogue Number
Mains adapter, 230V ac, UK plug	SM27/1
Mains adapter, 230Vac, 2 pin European plug	SM27/2
Adapter for car cigarette plug	SM27/3

For a comprehensive list of parts required by service engineers conducting internal repairs, please contact the sales department of Bibby Sterilin quoting both the model and serial number.

Repairs

Only spare parts supplied or specified by Bibby Sterilin or its agents should be used. Fitting of non-approved parts may affect the performance and safety features designed in to the instrument.

Any repairs or replacement of parts MUST be undertaken by suitably qualified personnel.

If in any doubt please contact the Technical Department at Bibby Sterilin at the address shown on the back cover, or the point of sale.

Bibby Sterilin Ltd.
Stone
Staffordshire
ST15 0SA
United Kingdom
Tel: +44 (0) 1785 812121
Fax: +44 (0) 1785 813748

Technical Specification

Max Speed	1300rpm
Capacity litres H ₂ O	1.5
Overall dimensions (w x d x h)	150 x 160 x 70mm
Net weight	1.3kg
Electricity supply	4 x alkaline batteries or 230V, 50 Hz
Battery type	1.5V, size D, ICE No. LR20
Battery life	6000 hours continuous use

Warranty

Bibby Sterilin Ltd. warrants this instrument to be free from defects in material and workmanship, when used under normal laboratory conditions, for a period of two (2) years. In the event of a justified claim Bibby Sterilin will replace any defective component or replace the unit free of charge.

This warranty does NOT apply if damage is caused by fire, accident, misuse, neglect, incorrect adjustment or repair, damage caused by incorrect installation, adaptation, modification, fitting of non-approved parts or repair by unauthorised personnel.

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These products meet the relevant EC harmonised standards for radio frequency interference and may be expected not to interfere with, or be affected by, other equipment with similar qualifications. We cannot be sure that other equipment used in their vicinity will meet these standards

and we cannot guarantee that interference will not occur in practice. Where there is a possibility that injury, damage or loss might occur if equipment malfunctions due to radio frequency interference, or for general advice before use, please contact the Technical Service Department of Bibby Sterilin Ltd.

Declaration of Conformity

Battery Operated Stirrer

Model SM27

This Stuart Scientific product complies with the requirements of the EU Directives listed below:

89/336/EEC Electromagnetic Compatibility Directive, amended by 92/31/EEC
and 93/68/EEC.

Compliance with the requirements of these Directives is claimed by meeting the following standards:

EN55014-1:1997 Electromagnetic compatibility Emission requirements for
household appliances electric tools and similar apparatus .
Product Family Standard.

EN55104:1995 Electromagnetic Compatibility. Immunity requirements for
household appliances, tools and similar apparatus. Product Family Standard.

Full Report Ref. No 19/98

The technical documentation required to demonstrate that the products meet the requirements of these Standards is available for inspection on request.

CE Mark Affixed 98

Signed:  (Mr C. R. Baggaley)

Date: 18/3/04

Authority: Technical Manager

for

Bibby Sterilin Ltd - Stone - Staffs - ST15 0SA - UK
Tel: +44 (0) 1785 812121 - Fax +44 (0) 1785 813748

INSPECTION REPORT

MODEL

SM4

SM27

ELECTRICAL SAFETY

1. Earth continuity
2. Insulation
3. Flash test

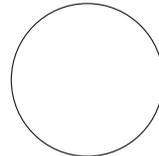
FUNCTIONAL

1. Indicators
2. Speed control
3. Visual acceptance

QUALITY CONTROL INSPECTOR

Bibby Sterilin

Bibby Sterilin Ltd

Beacon Road Stone
Staffordshire ST15 0SA
United Kingdom
Tel: +44 (0)1785 812121
Fax: +44 (0)1785 813748
www.bibby-sterilin.com

Bibby Sterilin France

BP 79-77793 Nemours Cedex
France
Tel: +33 1 64 45 13 13
Fax: +33 1 64 45 13 00
bibby@bibby-sterilin.fr

Bibby Sterilin Italia

Via Alcide de Gasperi 56
20077 Riozzo Di Cerro Al Lambro
Milano
Italy
Tel: +392 983 8969
Fax: +392 982 30211

Afora S.A.

Calle Aribau 240
08006 Barcelona
Spain
Tel: +343 93-306 98 00
Fax: +343 93-306 98 23
e-mail: marketing@afora.com
www.afora.com