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Use your scale with care, gently loading items to be weighed onto the center of the pan. Do not immerse the scale in water or other liquids, and avoid splashing the scale. Clean using a soft cloth dampened with water.

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Install 3-AA batteries (provided), or connect the AC adapter (optional accessory) to the power input jack at the rear of the scale. Do not apply excessive force or press on the pan.

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1

Weighing / Pesar / Pesage

Place sample on the pan and read the displayed result. The * indicator appears when the reading is stable. If using a container, press **ON/ZERO/Off** to tare the container before adding material.

If the applied load exceeds the capacity of the scale, an E will appear on the display and the load should be removed immediately.

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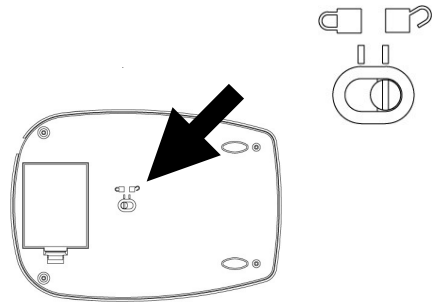
5

Slide the load cell lock to the unlock position for use. Lock for stacking, storing, or transport.

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2

Cumulative Weighing / Pesaje acumulativo / Pesage cumulatif

Place an empty container on the pan, allow the reading to stabilize then press **ON/ZERO/Off** to zero (tare) the weight of the container. The scale now displays zero and will display the weight of material added. Once the desired quantity is reached, press **ON/ZERO/Off** again to reset the display to zero before adding the next item.

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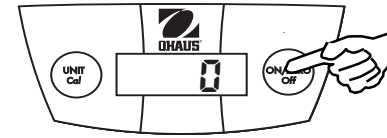
6

ON / ZERO / Off

Press this key to turn the scale on. This key is also used to zero the display. Press and hold this key to turn the scale off.

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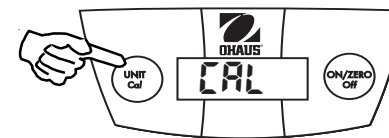
Refer to Optional Accessories for the correct calibration weights.

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
1. Press and hold **UNIT/Cal** until "CAL" is displayed, then release it.

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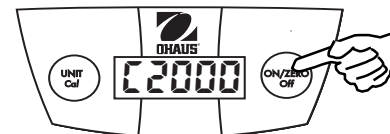2. Press **ON/ZERO**. The required weight will appear in grams.

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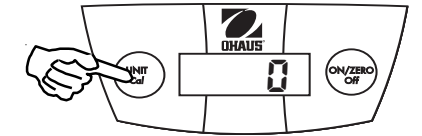
7

UNIT/Cal

Press this key to change the weighing unit. Press and hold to begin calibration.

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4

3. Place the appropriate weight on the pan, then press **ON/ZERO**.

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4. "PASS" will be displayed when calibration is complete.

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Note: If "CAL E" is displayed a calibration error has occurred (for example, an incorrect weight was used or the pan was disturbed). Repeat the procedure.

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8

Specifications

Model	CL5000F
Capacity x Readability	5000 x 1g 11lbs: 0.4 oz x 0.1 oz 176 x 0.1 oz
Repeatability (Std. Dev.) (g)	1
Linearity (g)	±1
Weighing Units	g, lb:oz, oz
Tare range	to capacity by subtraction
Stabilization time	2 seconds
Optimum operating temp.	65° - 77°F/18° - 25°C
Power requirements	3-AA alkaline batteries (included) AC adapter (optional accessory)
Battery Life	300 hours
Auto Shut-off	4 minutes no activity
Overload indication	⌈⌋

Optional Accessories	Ohaus Part No.
AC Adapter, 120V U.S.	90524-66
AC Adapter, 230V Euro	90524-63
AC Adapter, 240V U.K.	90524-64
AC Adapter, 240V Australia	72187149
Stainless steel pan cover	80010623
Carrying case	80010624
2000g Calibration mass	51026-16

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10

LIMITED WARRANTY

Ohaus products are warranted against defects in materials and workmanship for duration of the warranty. During the warranty period Ohaus will repair, or, at its option, replace any component(s) that proves to be defective at no charge, provided that the product is returned, freight prepaid, to Ohaus. This warranty does not apply if the product has been damaged by accident or misuse, exposed to radioactive or corrosive materials, has foreign material penetrating to the inside of the product, or as a result of service or modification by other than Ohaus. In lieu of product registration, the warranty period shall begin on the date of shipment to the authorized dealer. No other express or implied warranty is given by Ohaus Corporation. Ohaus Corporation shall not be liable for any consequential damages. As warranty legislation differs from state to state and country to country, please contact Ohaus or your local Ohaus dealer for further details.

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