# IKA

## IKA C 5010 / 5012

Betriebsanleitung	DE	3
Operating instructions	EN	16
Mode d'emploi	FR	29
Инструкция по эксплуатации	RU	42
使用说明	ZH	55

Indicaciones de seguridad	ES	68
Veiligheidsinstructies	NL	70
Norme di sicurezza	IT	73
S�akerhetsanvisningar	SV	75
Sikkerhedshenvisninger	DA	77
Sikkerhedsanvisninger	NO	80
Turvallisuusohjeet	FI	82
Normas de seguran�a	PT	84
�ποδείξεις ασφαλείας	EL	87
Wskaz�wki bezpiecze�stwa	PL	89
Bezpe�nostn� upozorn�n�	CS	91
Biztons�gi utasítások	HU	94
Varnostna navodila	SL	96
Bezpe�nostn� pokyny	SK	99
Ohutusjuhised	ET	101
Dro�ibas Nor�des	LV	103
Saugos Reikalavimai	LT	106
Инструкции за безопасност	BG	108
Indica�ii de siguran�a	RO	110

IKA - Werke, Germany  
Reg.-No. 004343

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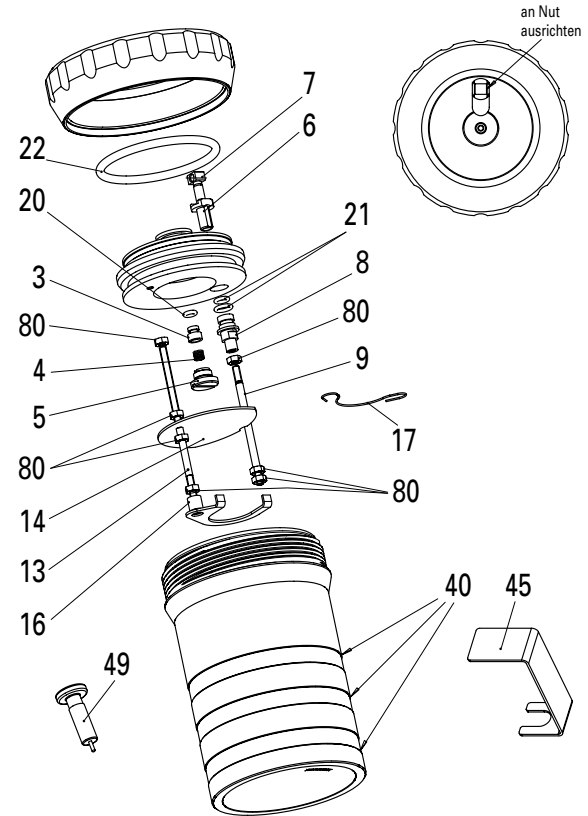
Source language: German

## Contents

	Page
CE Declaration of conformity	16
Warranty	16
Explanation of symbols	16
Safety instructions	17
Delivery scope C 5010/C 5012	19
Coding of the decomposition vessel	20
Opening the decomposition vessel	20
Fastening cotton thread	21
Closing the decomposition vessel	21
Combustion procedure	22
Cleaning the decomposition vessel	22
Halogens C 5012	22
Replacing the ignition wire	22
Care and maintenance	24
Replacing the crucible holder	24
Replacing the ignition electrode	24
C 5010 oxygen valve maintenance	24
C 5012 Oxygen valve maintenance	25
Replacing the C 5010 sealing ring	25
Leakage test	26
Degas	26
Decomposition vessel identificaton	26
C 5010 Technical data	27
C 5010 Spare parts list	27
C 5010 Spare parts diagram	27
C 5012 Technical data	28
C 5012 Spare parts list	28
C 5012 Spare parts diagram	28

EN

## CE Declaration of conformity

We declare under our sole responsibility that this product corresponds to the regulations for pressure vessels 2014/68/EU.

A copy of the complete EU Declaration of Conformity can be requested at [sales@ika.com](mailto:sales@ika.com).

## Warranty

In accordance with **IKA** warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our factory, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover worn out parts, nor does it apply to faults resulting from improper use, insufficient care or maintenance not carried out in accordance with the instructions in this operating manual.

## Explanation of symbols

This symbol identifies information **that is absolutely essential for your safety and health**. Failure to observe this information may result in negative effects for your health and injuries.

## Safety instructions

**Read the operating instructions in full before starting up and follow the safety instructions.**

- Keep the operating instructions in a place where they can be accessed by everyone.
- Ensure that only trained staff work with the appliance.
- Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.

### Intended purpose:

The C 5010 and C 5012 decomposition vessels must only be used for determining the gross calorific value of solid and liquid combustible materials in accordance with DIN 51900, BS 1016 T5, ISO 1928, ASTM 5468, ASTM 5865, and ASTM 4809.

Operation of decomposition vessels is only permitted in combination with the **IKA** calorimeter C 200, C 2000 or C 5000.

### Operating requirements:

The maximum **amount of energy input** into the decomposition vessel must not exceed **40000 J**. (Select the weight of the sample accordingly). The permissible **operating pressure** of **230 bar (23 MPa)** must not be exceeded. The maximum permissible **operating temperature** must not exceed **50 ° C**.

Do not fill the decomposition vessel too full of the sample. Only fill the decomposition vessel with oxygen up to a maximum pressure of **40 bar (4 MPa)**. Monitor the adjusted pressure on the pressure reducer.

Perform a check before every combustion to ensure there are no leaks (see the chapter entitled „Leakage test“).

### Explosive substances:

Many substances tend to combust in an explosive manner (for example because of the formation of peroxide). This may cause the decomposition vessel to burst.

**The C 5010 and C 5012 standard decomposition vessels must not be used for examinations on samples that are capable of exploding. It is absolutely essential to use a special high-pressure decomposition vessel to contain the sample in these cases!**

### Notes on the sample:

Substances of which the combustion behavior is not known must be examined for their combustion behavior before combustion in the decomposition vessel C 5010 and C5012 (danger of explosion).

If you are burning **unknown samples**, leave the room or **keep a safe distance** between you and the calorimeter

Benzoic acid must only be burned in the form of pellets!

Combustible dust and powder must be compressed into pellets before combustion. Oven-dry dust and powder such as wood chips, hay, straw, etc. burn in an explosive manner! They must be moistened first! Readily combustible liquids with a low vapor pressure must not be come in direct contact with the cotton thread (for example tetramethyl dihydrogen disiloxan)!

### Combustion residue, auxiliary materials:

In addition, toxic residues of combustion are possible in the form of gases, ash or precipitates on the inner wall of the decomposition vessel, for example

**Observe the accident prevention requirements applicable to the activity and the work station. Wear personal safety equipment.**

When handling combustion samples, combustion residues and auxiliary materials, the appropriate safety requirements must be observed. The following are examples of substances that may cause dangers: corrosive, easily flammable, capable of exploding, contaminated with bacteria, toxic.

#### **Oxygen:**

When working with oxygen, observe the appropriate requirements.

**Danger warning:** As a compressed gas, oxygen promotes combustion, supports combustion intensively and may react violently with combustible substances. **Do not use any oil or grease!**

#### **Using a crucible made of stainless steel:**

When using crucibles made of stainless steel, their condition should be carefully checked after every experiment.

A reduction in the thickness of the material may cause the crucible to burn and may damage the C 5010 and C5012 decomposition vessel.

For reasons of safety, crucibles must not be used any more after a maximum of 25 combustion procedures.

#### **Specification of the decomposition vessel:**

The decomposition vessel is manufactured in accordance with the regulation for pressure vessels 2014/68/EU. This can be recognized from the **CE symbol** with the identifying number of the testing station named. The decomposition vessel is a pressure device of Category III. The decomposition vessel has been subjected to an EC prototype test. The EU declaration of conformity represents our guarantee to you that this decomposition vessel complies with the pressure device described in the EC prototype test certificate. The decomposition vessel has been subjected to a pressure test at a **test pressure of 330 bar (33MPa)** and a leak test with oxygen at 30 bar (3MPa).

C5010 and C5012 decomposition vessels are **experiment autoclaves** and must be tested by a **professionally trained person** each time before they are used.

An individual application is understood here to mean a series of experiments that are performed under roughly the same conditions in terms of pressure and temperature. Experiment autoclaves must be operated in special chambers (IKA C 200, C 5000 and C 2000).

#### **Repeated tests:**

The decomposition vessel must be subject to repeated tests (internal tests and pressure tests) by a **person with professional training**. The intervals between tests must be determined by the operator based on experience, operating manner and the material used in the decomposition vessel.

**The declaration of conformity loses its validity if mechanical modifications are made to the experiment autoclaves or if stability can no longer be guaranteed as a result of heavy corrosion (for example holes eaten in it by halogens).**

The **threading** on the body of the decomposition vessel and cap screw in particular are subject to a high level of mechanical stress and must therefore be monitored regularly for **wear and tear**.

The condition of the seals must be checked for functionality must be ensured by means of a test for leaks (see the chapter entitled „Leakage test“)

If maintenance tasks, and especially the pressure testing, is not performed or is performed incorrectly, there is a risk the pressure vessel may burst or an uncontrolled internal fire may occur at the electrodes which could burn away the seals (oxyacetylene torch effect), thus posing a risk to life and limb.

Pressure tests and service tasks on the decomposition vessel must only be performed by **persons with professional training**.

**We recommend that the decomposition vessel be sent into our factory for inspection and repairs if necessary after either 1000 experiments or after one year or, depending on the application, even sooner than this.**

#### **Definition of person with professional training**

A person with professional training as defined in these operating instructions is someone:

1. whose training, knowledge and experience gained through practical activities ensures that that person will perform the tests in a proper manner.
2. who is sufficiently reliable.
3. who is not subject to any instructions in terms of testing activity.
4. who is equipped with suitable testing equipment if necessary.
5. who can provide suitable proof demonstrating compliance with the requirements listed in 1.

#### **Operating pressure containers:**

National regulations and laws for operating pressure containers must be observed!

Anyone who operates a pressure container must maintain it in proper condition, must monitor it and perform necessary maintenance and repair tasks without delay, and must take measures appropriate for the circumstances to ensure safety.

**A pressure container must not be operated if it exhibits defects that could endanger those working with it or third parties.**

You can obtain a copy of the pressure vessel regulation from Carl Heymann Press or Beuth Press.

**Use only original spare parts from IKA!**

## **Delivery scope C 5010/ C 5012**

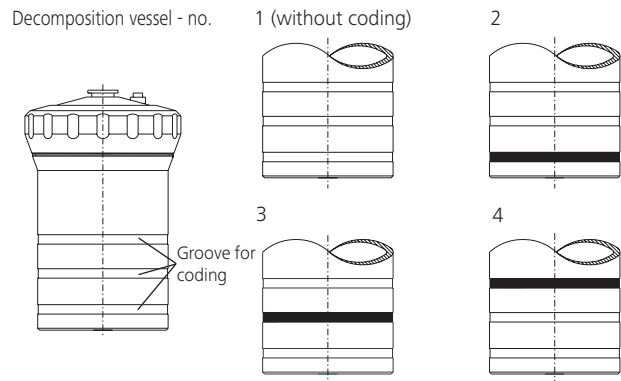
- 1 x decomposition vessel
- 1 x ignition wire
- 1 x O-Ring-Set
- 1 x C710.4 cotton thread
- 1 x gripp
- 1 x set of labels
- 1 x C 4 combustion crucible
- 2 x AF 5,5 double open ended wrench
- 1 x emergency-vent knob
- 1 x operating instructions

only C 5012:

- 1 x snap ring
- 1 x pressure spring


## Coding of the decomposition vessel

The label set included in the delivery quantity is used for coding. The individual strips are glued into the respective grooves around the decomposition vessel.


## Opening the decomposition vessel

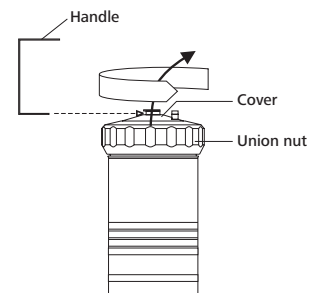- Unscrew union nut
- Use the handle in order to pull out the cover.


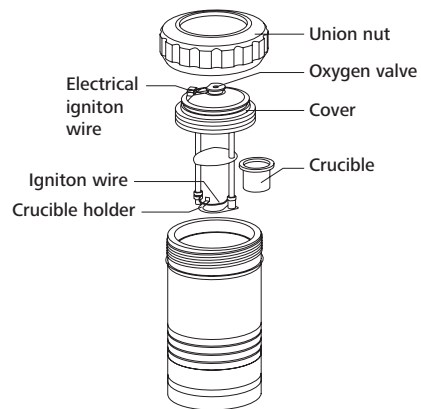
The decomposition vessel consists of the following individual parts:

## Fastening cotton thread

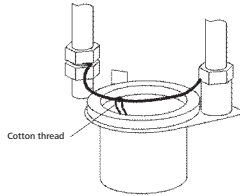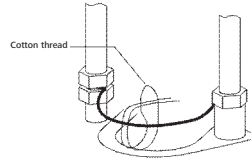
The decomposition vessel is equipped with a fixed ignition wire. A cotton thread has to be fastened to the ignition wire in order to carry out combustion.

Fasten cotton thread as shown in the diagram.  
Note: Beware the danger of burn-off in the case of incorrectly laid ignition wires!

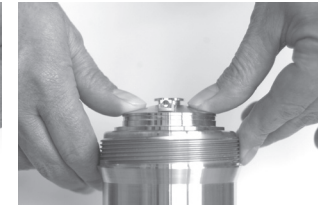
C 5010


C 5012

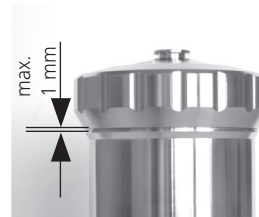- Place the union nut onto the lower section and tighten by hand.

C 5010

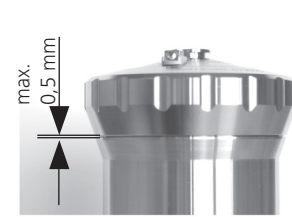C 5012

## Closing the decomposition vessel

- Place the cover onto the lower section and push down until it presses against the stop piece in the lower section.

C 5010

C 5012

- The decomposition vessel can now be filled.

## Combustion procedure

For the decomposition vessel charging procedure with fuel samples and the general procedure for performing combustion please refer to the relevant standards and the operating instructions for the **IKA C 5000 control / duo-control calorimeter system** or C 2000 and C 200.

When working with the **IKA C 5000 und C 2000 systems**, filling of the decomposition vessel takes place in the device itself; a filling station is used with the C 200 and C 248.

For the C 5000, venting of the decomposition vessel takes place in the device itself; for the C 2000 and C 200 the venting button **supplied with the device** is required.

## Cleaning the decomposition vessel

After each combustion the condensate has to be cleaned from all of the parts of the decomposition vessel which come into contact with the combustion products. It suffices to wipe out the parts with an absorbent, non-fibrous cloth. Combustion residues in the crucible, e.g. carbon black or ash are removed in the same manner.

The decomposition vessel may not be mechanically cleaned under any circumstances if it is not capable of being cleaned in the manner described (e.g. due to burn-in or pitting).

In such cases please contact your **competent service center** or send in the decomposition vessel to the factory for inspection and cleaning.

## Halogens C 5012

Pitting which severely impairs operator safety may be caused on the inner surfaces of the decomposition vessel from combustion products which contain halogen. If the halogen content of the samples is significantly greater than the trace range, then decomposition vessel C 5012, which is specially designed for such applications, shall be used.

To enable the **quantitative evaluation of halogens and sulphur**, the C 5012 decomposition vessel is preconditioned before shipment. This means that the inner surfaces and the fittings of the decomposition vessel are catalytically activated by means of a patented procedure..

Parts thus treated are recognizable because of their opaque, tarnished and partly spotty surfaces. This appearance of this surface is intentional and should not be confused with traces of wear.

This state is important for the following analysis and may not be treated with brushes or other hard materials. Wiping and drying should be done using a soft cloth.

**The C 5010 decomposition vessel may not be used for such applications because of the safety risk and insufficient analysis results.**

## Replacing the ignition wire

The C 5010 and C 5012 decomposition vessels are equipped with a fixed ignition wire. Since the ignition wire is subjected to a high degree of material stress during combustion, it should be replaced after about 100 tests at the latest. However, this largely depends on the type of samples fired.

The ignition wire must also be replaced if corrosion occurs, in order to prevent drift in the results.

The ignition wire from the C 5010 and C 5012 decomposition vessels are not interchangeable because of the difference in the alloys.

Leakages of the decomposition vessel are the results (see the sections entitled „Leadkage test“).

**ATTENTION!**

**Leakages at the decomposition vessel container can cause heavy damage to the calorimeter and the decomposition vessel. Ignoring can lead to health impairment and injury.**

#### **Detaching the ignition wire on the ignition electrode (item 20):**

- Fix the ignition electrode (item 20) in place at the spanner surfaces using the AF 5.5 double open-ended spanner supplied.
- This will prevent the ignition electrode from being twisted.
- Loosen the nut (item 30) using the AF5 .5 double open-ended spanner.

#### **Detaching the ignition wire on the crucible holder (item 25):**

- Firmly hold the crucible holder and deflector and unscrew the hexagonal nut using the AF 5,5 double open-ended wrench.

**Note: The crucible holder should not be twisted while doing so.**

#### **Insert / Tightening of the ignition wire:**

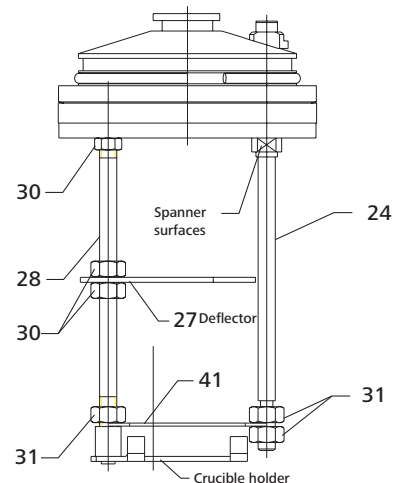
The ignition wire is shaped in such a way that the electrodes of the decomposition vessel fit into the side holes.

The following points have to be strictly observed:

- The ignition wire has already been shaped in such a way that that it lies above the crucible holder (note curvature).
- Align the ignition wire horizontally between the electrodes.
- Do not press the electrodes together. The deflector of the crucible holder may be come into contact with the ignition electrode.
- The ignition electrode must not be twisted and thereby loosened.

- Follow the sequence in reverse to tighten the nut.
- Ensure that there is a sound contact with the ignition wire.

Note: Beware the danger of burn-off in the case of incorrectly laid ignition wires!

Electrode with ignition wire

#### **C 5010**

- Pos. 24 Electrical ignition wire
- Pos. 27 Deflector
- Pos. 28 Shaft
- Pos. 30 Hexagonal nut M4 A2
- Pos. 31 Hexagonal nut M3 Hastelloy
- Pos. 41 Ignition wire

#### **C 5012**

- Pos. 9 Electrical ignition wire
- Pos. 13 Shaft
- Pos. 14 Deflector
- Pos. 17 Ignition wire
- Pos. 80 Hexagonal nut M3 Hastelloy

**ATTENTION!**

**The use of alternative materials can pose a risk to life and limb in the case of the Hastelloy nuts (Pos. 30).**

## Care and maintenance

The work described in the following may only be performed by personnel with professional training or **IKA**.

All work on the fittings of the decomposition vessel require a leak test and new pressure testing.

Furthermore, reference is made to the regulation for pressure vessels 2014/68/EU.

Send in the case of service the equipment back in the original packaging. Storage packing is not sufficient for the back dispatch. Use additionally a suitable transportation packing.

## Replacing the crucible holder

- Remove ignition wire
- Firmly hold crucible holder and unscrew the hexagonal nut at the upper end of the electrode in counterclockwise direction.
- Screw the crucible holder out of the cover.
- Insert new crucible holder and align it in such a way that the slot in the deflector is flush with the ignition electrode located opposite (crucible holder toward the inside).
- Tighten locknut.

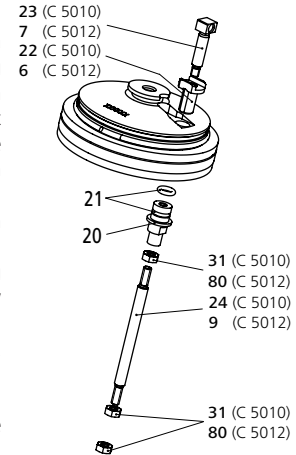
## Replacing the ignition electrode

- Remove ignition wire.
- Unscrew electrode (item 20) in a counterclockwise direction using the double open-ended wrench and holding up the female insert contact (item 23) using a suitable tool (item 20 screwed together with item 23).
- Remove individual parts as shown in the following diagram.
- A new O-ring (item 21) (cf. O-ring set) should be used when the new electrode is inserted.
- Join individual parts (items 20, 22 and 23) and screw together firmly.

**Note!**

After being tightened the female insert contact (item 23) and insulating bushing (item 22) have to be flush with the recess in the cover. Repeat procedure if required.

## C 5010 oxygen valve maintenance

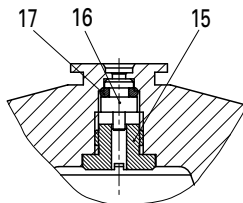
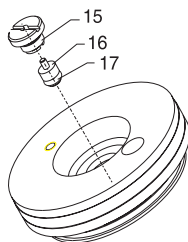
The oxygen valve is used to fill and empty the decomposition vessel. The O-ring sealing may wear because of aggressive combustion gases, which means that the decomposition vessel may not be tight.

In order to replace the O-ring the valve inside the cover has to be opened:

- Unscrew locking screw (item 15) using a slotted screwdriver.
- Remove valve piston (item 16) and O-ring (item 17).
- Replace O-ring (cf. O-ring set).
- Insert piston in such a way that the guide pin point toward the locking screw.
- Tighten locking screw.

Mounted oxygen valve:


## C 5012 oxygen valve maintenance

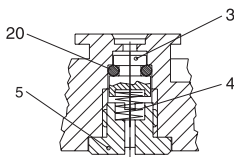Disassembly of the oxygen valve for the and C 5012 decomposition vessel is identical to the procedure for the C 5010 decomposition vessel. However, the components differ in their design.

- Pos. 3 valve piston
- Pos. 4 pressure spring
- Pos. 5 locking screw
- Pos. 20 O-Ring

Assembly of the components shall be performed as specified in the preceding diagram.

The pressure spring (item 4) does thus the oxygen valve the decomposition vessel is self-sealing.

## Replacing C 5010 sealing ring

In order to raise the sealing ring (item 13) the circlip (item 11) has to first be pried out of the groove in the cover.

This is performed using a medium-sized slotted screwdriver, the flat side of which is pressed between the circlip and the cover. When the screwdriver is properly positioned the circlip is removed by pushing back the cover edge.

Now the compression ring (item 12) and the sealing ring (item 13) may be lifted.

Pos. 11 circlip

Pos. 12 compression ring

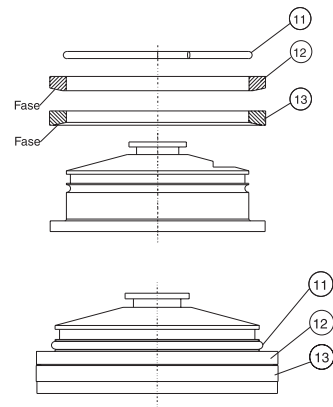
Pos. 13 sealing ring

When joining the individual parts make certain that the chamfers of the compression ring and sealing ring point toward the side of the cover (see diagram).

Mounted components:

The circlip (item 11) has to snap exactly into the cover groove. There is hardly any play between the circlip and the compression ring after assembly.

The circlip may not be put into the upper groove of the cover groove because the decomposition vessel would not be tight.

A leak test and pressure testing shall be performed after assembly has been successfully completed (see the sections entitled „Leakage test“).

## Leakage test

**Perform a leakage test before every combustion process. Always check the lid is properly sealed with the decomposition vessel after carrying out maintenance. Follow the safety instructions.**

- Fill the decomposition chamber in the calorimeter system or using the C 248 oxygen filling station.
- Remove the decomposition vessel and immerse fully in a container filled with water.
- Check the decomposition vessel for oxygen leaks by watching for bubbles.
- If bubbles are seen then the decomposition vessel is not sealed correctly.
- Vent the decomposition vessel and check the seals.

**Never use a leaking decomposition vessel in the calorimeter!**

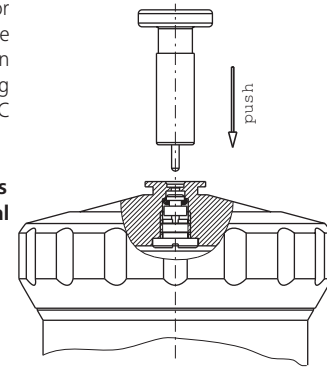
## Degase

For the C 5000, venting of the decomposition vessel takes place in the device itself; for the C 2000 and C 200 the venting button supplied with the device is required.

The vent knob is supplied for manual release of tension in the decomposition vessel. (Even in case of error function of the filling unit in the calorimeter C 2000 / C 5000).


**Escaping decomposition gases might be detrimental to health.**

## Decomposition vessel identification

The individual parts of the decomposition vessel are clearly identified by means of the engraved serial number and the technical data.

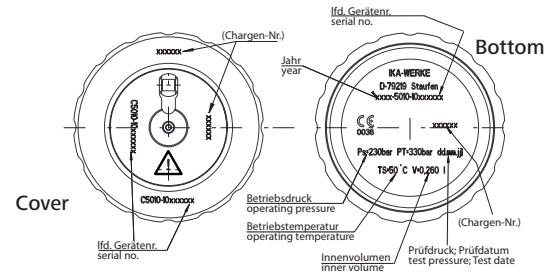The individual parts of several decomposition vessels (cover, union nut and bottom part) are not interchangeable.

If one of the parts indicated becomes damaged then the complete decomposition vessel has to be sent to the factory for testing.


## C 5010 Technical data

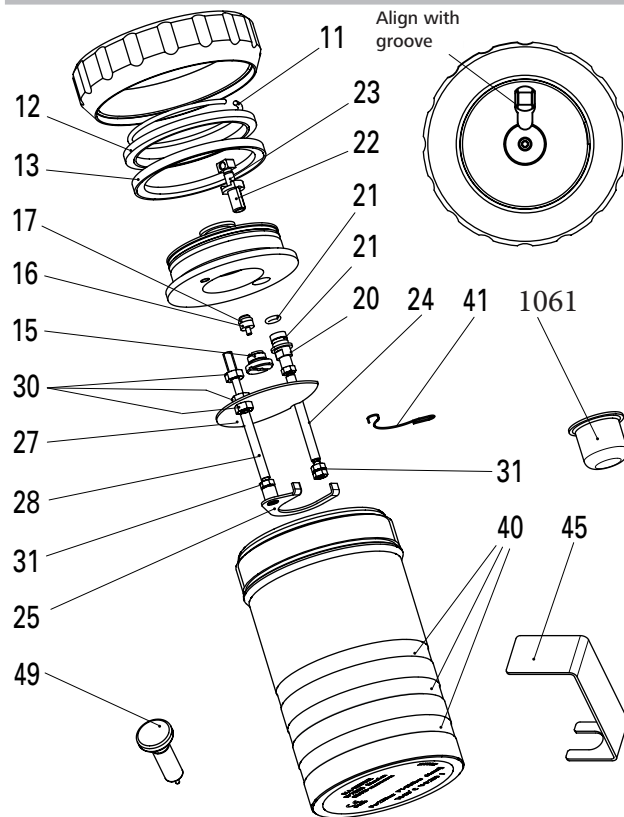Operating pressure	max. 230 bar (23 MPa)
Test pressure	330 bar (33 MPa)
Volume	260 ml
Nuclear temperatures	max. 1300 °C
Sample decomposition	isochor
Oxygen valve	plate valve
(vacuum operation not possible)	
Dish holder	as in DIN 51 900
Material	1.4439 2.4610.40

## C 5010 Spare parts list

Pos. 11 Spring ring	Pos. 25 Crucible holder
Pos. 12 Compression ring	Pos. 27 Deflector
Pos. 13 Sealing ring	Pos. 28 Shaft C 5010
Pos. 15 Locking screw	Pos. 30 Hexagonal nut M4 A2
Pos. 16 Valve piston	Pos. 31 Hexagonal nut M3 Hastelloy
Pos. 17 O-ring 3.1 x 1.6	Pos. 40 Lable set
Pos. 20 Bushing	Pos. 41 Ignition wire C5010.3
Pos. 21 O-ring 4 x 1.5	Pos. 1061 C4 quartz crucible
Pos. 22 Insulating bush	Pos. 45 Grip
Pos. 23 Female insert contact	Pos. 49 emergency-vent knob
Pos. 24 Ignition electrode	

## C 5010 Spare parts diagram

## C 5012 Technical data

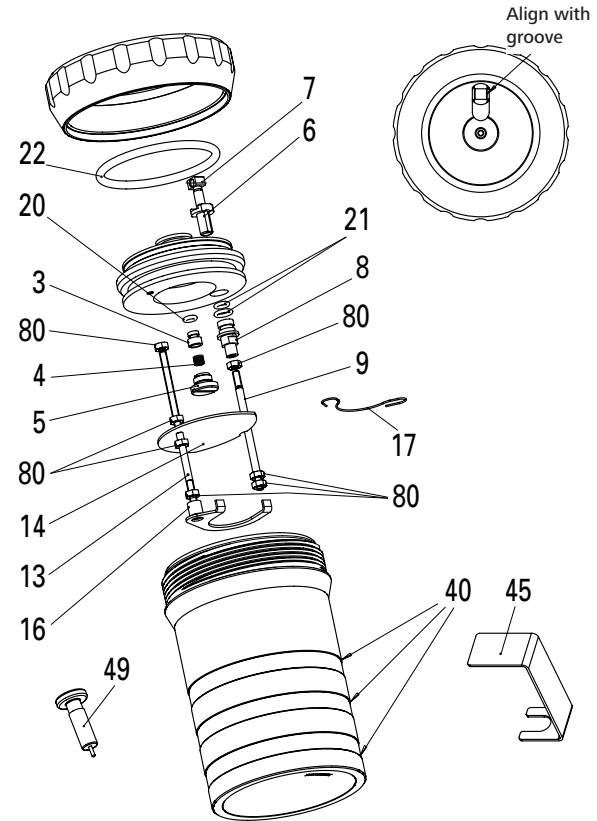
Operating pressure	max. 230 bar (23 MPa)
Test pressure	330 bar (33 MPa)
Volume	210 ml
Nuclear temperatures	max. 1300 °C
Sample decomposition	isochor
Oxygen valve	plate valve
(vacuum operation not possible)	
Dish holder	as in DIN 51 900
Material	2.4610.40 conditioned

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IKA Korea Ltd.
IKA India Private Limited
IKA England LTD.