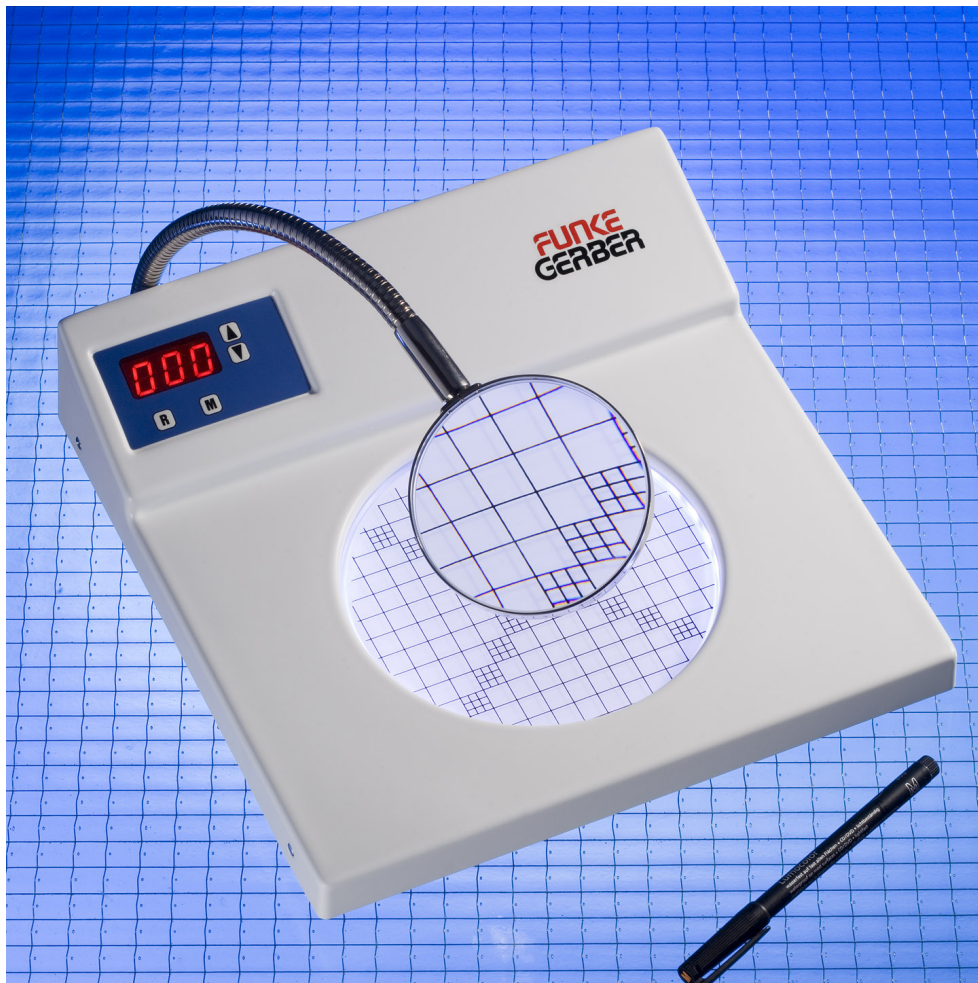
**FUNKE
GERBER**

ColonyStar

Art. No. 8500 - Deutsch

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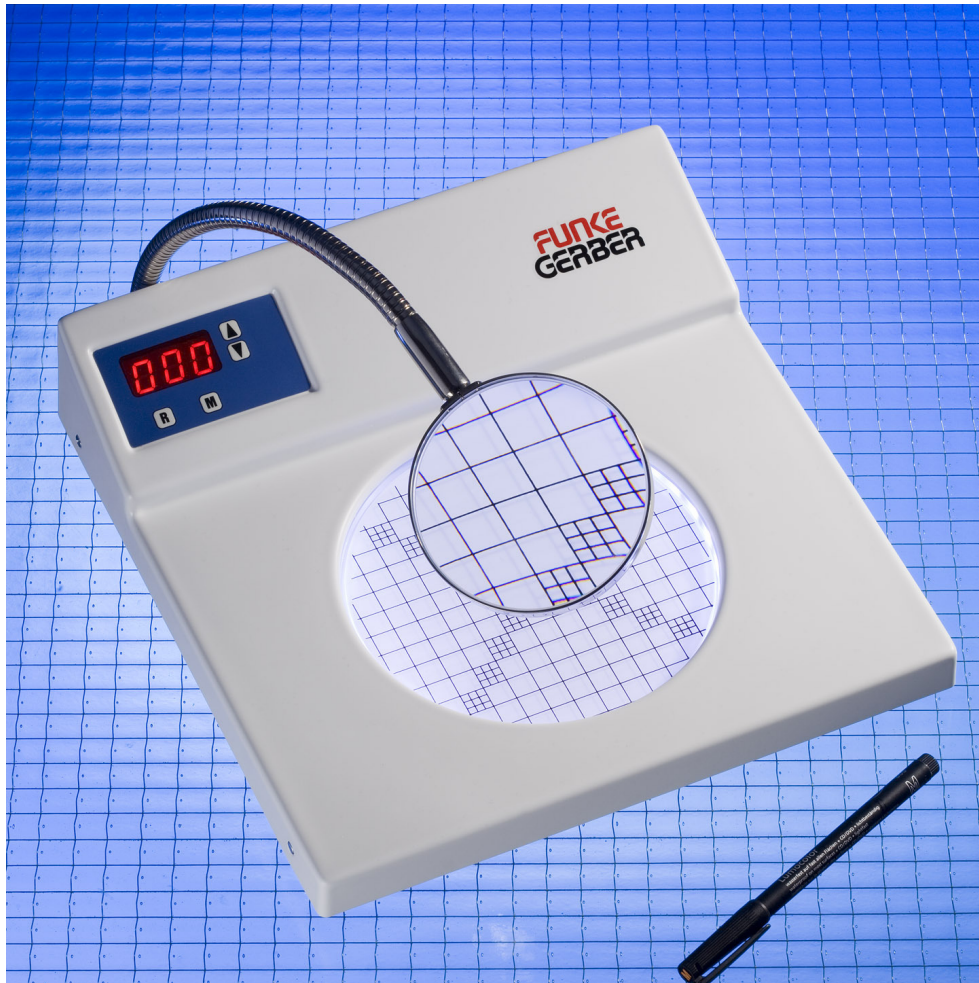
Safety Instructions

Please read and observe!

Please check if the voltage matches the details on the rating plate before connecting the unit to the mains.

The product should not be opened because of risk of electric shock. If the device must still be open in exceptional cases, pull out the power plug before opening the device

Put no heavy objects on the illuminated area or drop them on the area, this results in damage to the device.

Cleaning Instructions

The housing can be cleaned with a damp cloth. Care must be taken that no liquid enters the unit.

The illuminated area must not be cleaned with organic solvents, such as Thinners, alcohols, etc., this can lead to damage of the illuminated area.

Manual

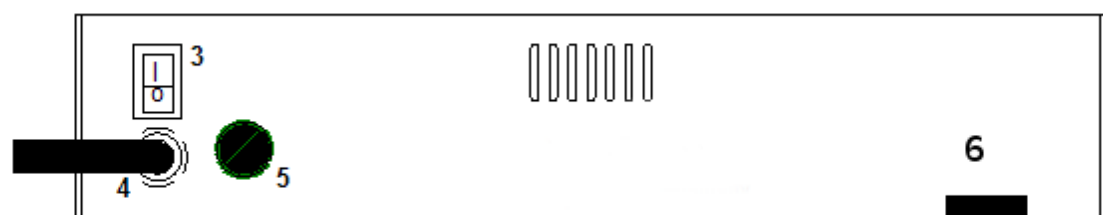
The device is used for easy and fast enumeration of Petri dishes. With the help of the pressure-sensitive light area and a marker pen it is easier to count substantially - by marking the respective colony with fiber pen, the electronic counter is increased by one.

Small Petri dishes are adapted with the help of the reducing insert.

With the frosted white milk glass plate the Petri dishes under direct lighting (illumination) are analyzed. An indirect, side lighting, important in the counting of Petri dishes containing medium light cultures, is achieved using the clear glass plate with dark field.

Delivery

- ColonyStar
- Marker pen
- Magnifying glass
- Frosted glass, white
- Clear glass with black side
- Reducing insert
- Manual

Device structure

- 1 counter and control unit
- 2 The glass can be replaced or removed by gently turning the device over.
- 3 Power switch to turn on and off the device
- 4 mains - see safety instructions
- 5 Fuse
- 6 Hole for the magnifying glass

The counting and control unit

Button R	reset button
Button M	Menu: selections of functions
Button ▲	Increase button
Button ▼	Decrease button

By repeated pressing the key **M** the modes are selected in the following order: **brightness** → **colour** → **sensitivity** → **signal generator** → **counting**

In counter mode, the counter can be reset to zero by pressing R. With ▲ and ▼ a manual correction of the displayed count is possible.

Mode brightness

By pressing ▲ and ▼ the brightness can be increased and decreased.

Mode colour

By pressing ▲ ▼ and the color can be changed.

Mode sensitivity:

From steps 0 (lowest sensitivity) to 9 (highest sensitivity) the sensitivity can be chosen by the keys ▲ and ▼

Modus signal generator:

The acoustic signal generator can be turned off and on by the ▲ and ▼ keys.

Setup values remain stored even after turning off the device and are restored when resuming operation.

The counter value is not stored.

Fault diagnosis

The illuminated area remains dark, even though the device was connected to the mains and switched on.

If the counts of the control unit lights, check if brightness is set in a mode value greater than zero. If the counting unit is dark too, check the fuse and replace it with a similar model.

The counter does not count, or only rarely when marking the colonies.

When the counter is not responding even at very high sensitivity to pressure, there is probably a defect of the device.

Technical data

Line voltage:	230 V AC / 50 Hz
Weight:	2.2 Kg
Height:	80 mm
Width:	275 mm
Depth:	300 mm
Maximum count:	999
Brightness levels:	65
Number of Colors:	65

Ordering

Description	Article
ColonyStar	8500
Frosted glass, white	8502-001
Clear glass with dark side	8505
Magnifying glass	8501-001
Reducer	8502-002
Marker pen	8504

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Z.A de Gesvrine – 4 rue Képler – B.P.4125
44241 La Chapelle-sur-Erdre Cedex – France
t. : +33 (0)2 40 93 53 53 | f. : +33 (0)2 40 93 41 00
commercial@humeau.com

EC Declaration of Conformity
under the EC Machine Directive 06/42/EG

We declare that the machine referred to herein below, as put on the market, has been designed and manufactured to the fundamental safety and health standards as defined in the EC Machine Directive. This Declaration shall become null and void in the case of any modification made without prior consent.


Designation of the machine:	Colony Star
machine typ:	colony counter
Serial number :	8500-
Applicable EC Directives:	EC Machine Directive 06/42/EC Low Voltage 2006/95/EC Electromagnetic Compatibility 2004/108/EC

Berlin, this
Funke-Dr.N.Gerber Labortechnik GmbH

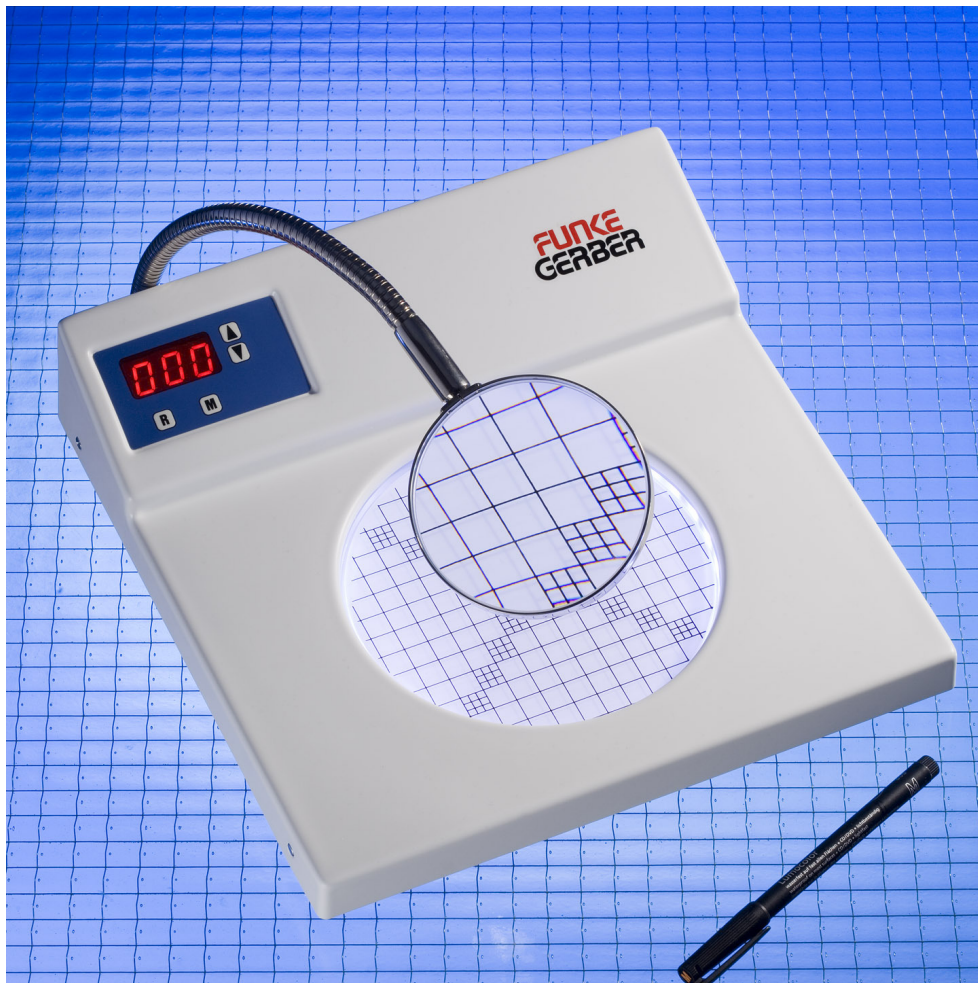Dipl.-Ing. K. Schäfer, Managing Director

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