
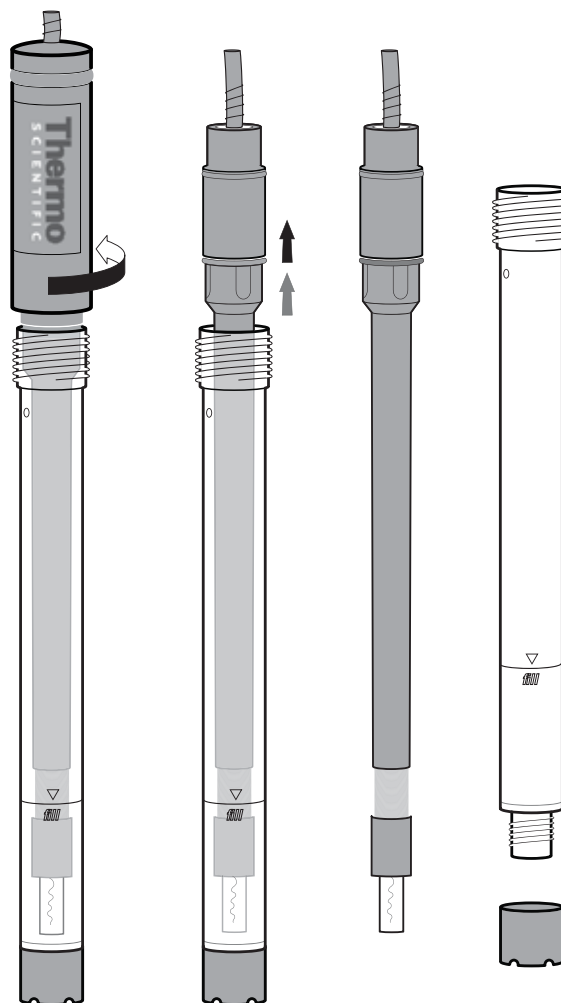
Thermo Scientific
Orion High Performance
Ammonia Electrode



Quick Start Guide

Guía rápida / Kurzanleitung / Guide de démarrage / Guida rapida

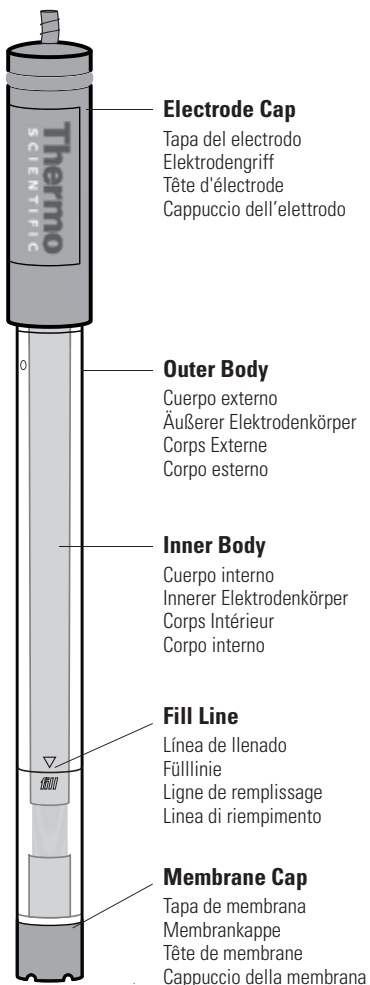Loose Membranes
(Cat. No. 951214)

Membranas sueltas
Lose Membranen
Membranes souples
Membrane di ricambio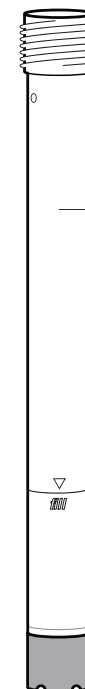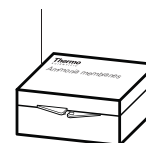

Preassembled Outer Body and Membrane Cap, includes outer body, membrane cap and membrane.
(Cat. No. 951215)

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Environmental Instruments
Water Analysis Instruments

166 Cummings Center
Beverly, MA 01915 USA

Toll Free: 1-800-225-1480
Tel: 1-978-232-6000
Dom. Fax: 1-978-232-6015
Int'l Fax: 978-232-6031
www.thermo.com/water

©2007 Thermo Fisher Scientific Inc.
All rights reserved.

254004-001 Rev.A 1007

Electrode Preparation

1. Unscrew the electrode cap from the electrode body. **See A.**
2. Remove the inner body from the outer body assembly and save the outer body assembly for spare parts. **See B.**
3. Soak the glass tip of the inner body in electrode filling solution (Cat. No. 951209) for at least two hours. **See C.**
4. Remove the preassembled outer body and membrane cap (Cat. No. 951215) from the electrode box. Verify that the membrane is not wrinkled or torn. **See D.**
5. Fill the preassembled outer body and membrane cap to the fill line with electrode filling solution. **See E.**
6. Insert the inner body into the preassembled outer body and membrane cap. Ensure that the inner body is fully inserted in the top of the outer body. **See F.**
7. Screw on the electrode cap until finger tight.
8. Rinse the electrode by immersing it in distilled water and blot it dry.
9. Soak the electrode in 100 mL of 1 ppm standard with 1 mL of alkaline reagent (Cat. No. 951011) for at least 15 minutes before use.

Electrode Storage

Between measurements— store the electrode in 100 mL of 1 ppm standard with 1 mL of alkaline reagent (Cat. No. 951011).

For overnight storage and storage up to one week— store the electrode in the electrode filling solution (Cat. No. 951209).

For storage over one week— disassemble the electrode and rinse all the components with distilled water. Dry and reassemble the electrode without electrode filling solution.

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