BEDIENUNGSANLEITUNG
OPERATING MANUAL
MODE D'EMPLOI
INSTRUCCIONES DE OPERACIÓN

ba15356defs03 03/2015

CN 500 DIN, CN 500 BNC
CN 800 DIN, CN 800 BNC

a xylem brand

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a xylem brand

**Note**

The latest version of the present operating manual can be found on the Internet under www.WTW.com. It is also available in other languages.

Copyright


© 2014, WTW GmbH

Reprinting - even as excerpts - is only allowed with the explicit written authorization of WTW GmbH.

Printed in Germany.

Overview

View

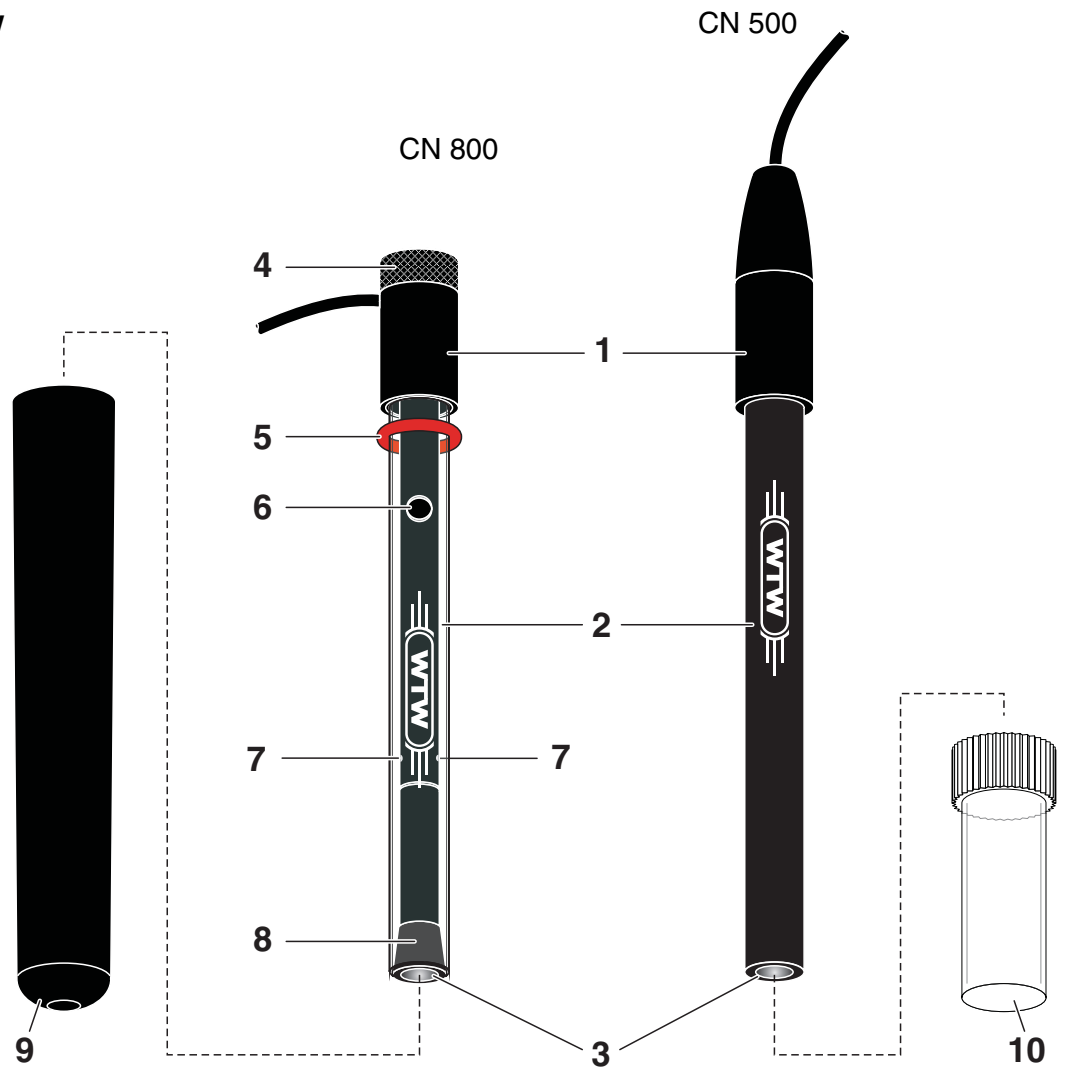1	Connection head with connection cable
2	Shaft
3	Membrane
4	Knurled nut
5	O-ring
6	Filling opening for bridge electrolyte
7	Inner junctions
8	Ground junction
9	Storage vessel for CN 800
10	Protective cap for CN 500

Commissioning

Danger

When dealing with acidic solutions that contain cyanide, there is danger of poisoning by released hydrocyanic acid. It is imperative to add the alkaline reagent MZ/NH₃/CN to any sample in order to raise the pH value of the sample to > 10 and prevent the release of hydrocyanic acid. Carry out cyanide determinations in acidic media with special safety measures only, e.g. in a fume hood.

Combination electrode CN 800

1	Remove the storage vessel.
2	Fill the bridge electrolyte ELY/BR/503 into the filling opening.
3	Shortly press the shaft of the combination electrode against the connection head to wet the ground junction with bridge electrolyte.
4	When doing so refill any spilled bridge electrolyte so that the inner junctions are covered with bridge electrolyte.
5	Rinse the combination electrode with deionized water.
6	Wipe the shaft using a clean paper towel.


Double rod electrode CN 500 + R 503

For measurements with the CN 500 electrode, a reference electrode is required (e.g. R 503). The two electrodes together form a double rod combination electrode.

1	Put the reference electrode into operation (see operating manual of the reference electrode).
2	Remove the protection cap of the electrode.
3	Rinse the electrode with deionized water.
4	Wipe the shaft using a clean paper towel.

Conditioning, calibration, measurement

General information

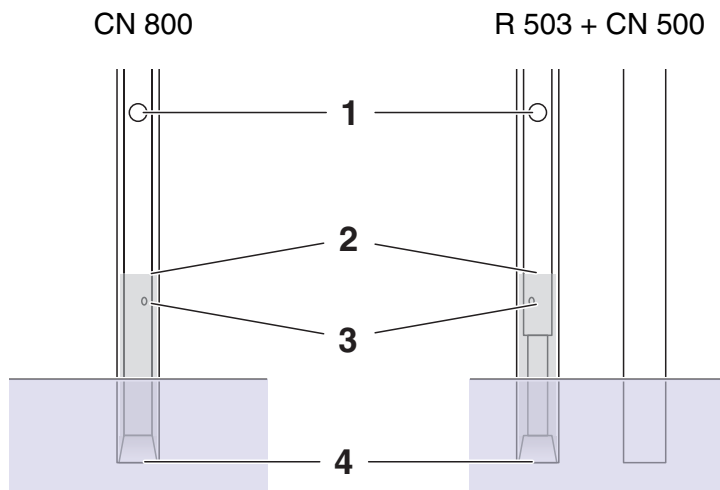When operating the electrode ensure that

- the filling opening (1) for the bridge electrolyte is open
- the inner junctions (3) are covered with bridge electrolyte
- no air bubbles are in the bridge electrolyte
- the depth of immersion is within the optimum range:

Minimum depth of immersion

The ground junction (4) must be covered

Maximum depth of immersion

Approx. 1 cm below the fluid level (2) of the bridge electrolyte

Before measuring

1	Condition every ion sensitive electrode in standard solution before using it. To achieve a good responding behavior, the concentration of the standard solution should be similar to the concentration expected of the test sample.
2	Remove any air bubbles in the bridge electrolyte by slightly knocking against the shaft.
3	Calibrate according to the operating manual of the meter and the analysis specification.

The conditioning time depends on the previous history of the ion sensitive electrode. Connect the electrode(s) to the meter and switch to the measured value display (see operating manual of the meter). The ion sensitive electrode is sufficiently conditioned when the measured value no longer changes considerably under otherwise constant conditions (temperature, composition of the solution).

NOTE

Avoid measuring samples with cyanide concentrations of more than

25 mg/l, as otherwise the electrode could be damaged irreversibly. If you expect the concentration to be higher, dilute the sample accordingly.

In the case of an erroneous calibration immediately after commissioning, polish the membrane surface (see MAINTENANCE).

Malfunctions can possibly occur due to the use of a magnetic stirrer. If the values are fluctuating heavily, we recommend to switch off the magnetic stirrer during measurement.

Sample preparation

Add 1 % MZ/NH₃/CN solution. This sample conditioning solution creates optimum conditions for measuring. It provides a constant ionic strength and similar diffusion potentials at the reference electrode in standard solution and test sample.

If you would like to have more detailed information concerning sample preparation and measuring procedures, WTW provides a large number of application reports for various applications.

Response times

The response time depends on the concentration range. It is

- several seconds at high concentrations,
- several minutes near the detection limit.

The measured value is stable if the value does not change by more than 0.1 mV within 30 seconds.

Interferences

- Complexes with Ni²⁺, Cd²⁺, Cu²⁺
- Strong reducing agents
- Interfering ions: 10 % error with the following concentration ratio (concentration ratio = interfering ion / measured ion):

Cl ⁻	I ⁻	Br ⁻	S ²⁻
1x10 ⁶	0.1	5x10 ³	must not be present

Aging

Please note that every electrode undergoes a natural aging process. The response time increases and the slope decreases with the age of the electrode. The following factors shorten the lifetime considerably:

- Unsuitable storing conditions (see section STORAGE)
- Special measuring conditions (e.g. organic solutions, frequent measurement with high concentrations of interfering ions)
- High temperatures
- High changes in temperature

The warranty does not cover damage caused by measuring conditions and mechanical damage.

Maintenance

- Combination electrodes:
Refill any used up bridge electrolyte.
- In the case of increased response times of the electrode perform the following activities:
 - Polish the membrane surface using an SF/K polishing strip (see section RECOMMENDED ACCESSORIES). To do so, moisten a polishing strip with deionized water, and with circular movements polish the membrane surface for approx. 30 seconds.
 - Then recondition the electrode (see section CONDITIONING, CALIBRATION, MEASUREMENT).
- Clean the inside of the combination electrode. To do so, open the combination electrode as follows:

Opening the combination electrode

1	Unscrew the knurled nut from the connection head.
2	Position the connection cable upward in a straight line with the connection head.
3	Unscrew the connection head.
4	Push the connection head and pressure spring over the connection cable.
5	Push the inner body out of the shaft.

NOTE

**Never pull the connection cable of the electrode.
The cable might be damaged.**

Storage

Between two measurements

Condition the electrode in standard solution.

Overnight

CN 500: Rinse the electrode with deionized water and store it in a dry place.

CN 800: Rinse the combination electrode with deionized water and dry it. Then put on the storage vessel and store the electrode upright.

Longer than overnight

CN 500: Rinse the electrode with deionized water, then dab it dry with a clean paper towel and put on the protective cap. Store the electrode in a dry place.

CN 800: Remove the bridge electrolyte, then rinse the combination electrode with deionized water and dab it dry with a clean paper towel. Store the combination electrode in the storage vessel in a dry place. For the following measurement, prepare the combination electrode as described in section COMMISSIONING.

Store the reference electrode according to the instructions in its operating manual.


Recommended accessories

Description	Model	Order no.
Reference electrode for electrode CN 500	R 503/P* R 503/D**	106570 106571
Bridge electrolyte	ELY/BR/503	106 575
ISA sample conditioning solution for CN ⁻ measurement	MZ/NH3/CN	150130
Polishing strips (24 pieces)	S/FK	180130

* Pin plug

** Banana plug

Calibration line of a cyanide sensitive electrode

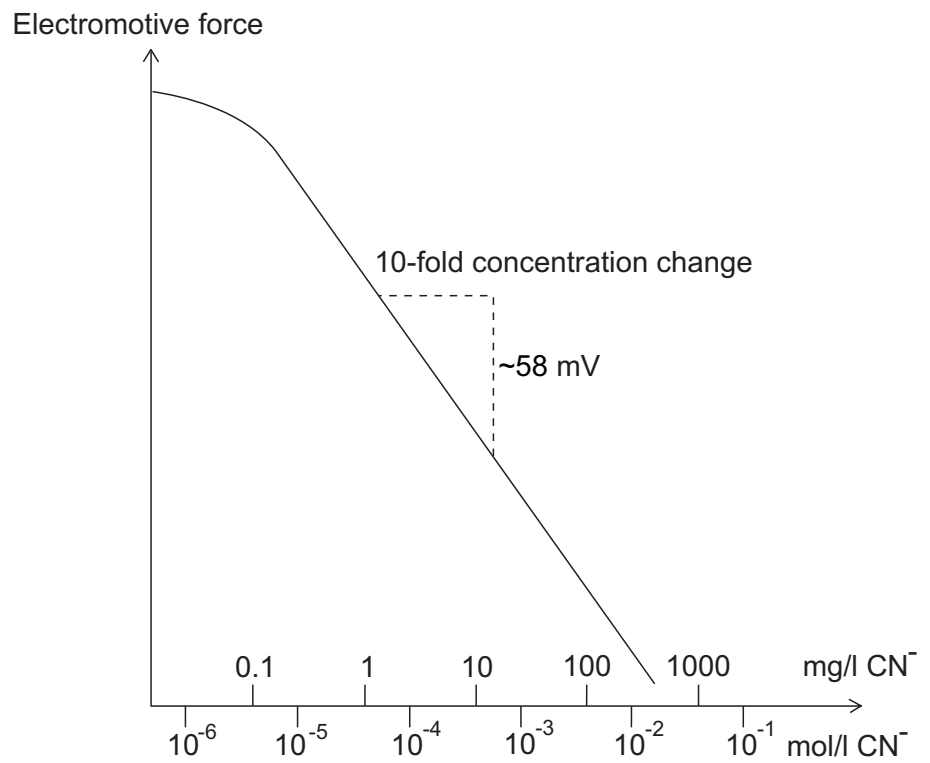What to do if ...

Measured value unstable	Cause	Remedy
	– Inner junctions not sufficiently wetted with bridge electrolyte (CN 800)	– Fill up bridge electrolyte until the inner junctions are covered with bridge electrolyte
– Inner junctions encrusted (CN 800)	– Leave the bridge electrolyte to react on the inner junctions for some hours until the crusts have dissolved.	
– Ground junction contaminated (CN 800)	– Rinse ground junction with bridge electrolyte	
– Cable broken	– Exchange the electrode	
Slope too low	Cause	Remedy
	– Membrane surface contaminated	– Polish the membrane surface with polishing strip S/FK (see MAINTENANCE)
– Conditioning time too short	– Extend the conditioning time	
– Standard solutions too old	– Use new standard solutions	
– Inner junctions encrusted (CN 800)	– Leave the bridge electrolyte to react on the inner junctions for some hours until the crusts have dissolved.	
– Electrode defective	– Exchange the electrode	

Technical data

Measuring ranges	0.2 ... 260 mg/l CN ⁻ (recommended 0.2 ... 25 mg/l CN ⁻) 8 x 10 ⁻⁶ ... 1 x 10 ⁻² mol/l CN ⁻ (recommended 8 x 10 ⁻⁶ ... 1 x 10 ⁻³ mol/l CN ⁻)	
Reproducibility	± 2 %	
pH range	0 ... 14	(Before measuring, adjust the pH value of the sample to pH 10 to 14)
Temperature range	0 ... 80 °C (shortly up to 100 °C)	
Resistance	< 30 MΩ	
Length	CN 500:	170 mm (including 50 mm connection head)
	CN 800:	153 mm (including 33 mm connection head)
Diameter	Shaft:	12 mm
	Connection head:	16 mm
Cable length	1 m	
Plug	DIN plug or BNC plug, depending on the design.	

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What can Xylem do for you?

We're a global team unified in a common purpose: creating innovative solutions to meet our world's water needs. Developing new technologies that will improve the way water is used, conserved, and re-used in the future is central to our work. We move, treat, analyze, and return water to the environment, and we help people use water efficiently, in their homes, buildings, factories and farms. In more than 150 countries, we have strong, long-standing relationships with customers who know us for our powerful combination of leading product brands and applications expertise, backed by a legacy of innovation.

For more information on how Xylem can help you, go to xylem.com.

Service address:

Xylem Analytics Germany
Sales GmbH & Co. KG
WTW
Dr.-Karl-Slevogt-Str. 1
82362 Weilheim
Germany

Tel.: +49 881 183-325
Fax: +49 881 183-414
E-Mail wtw.rma@xylem.com
Internet: www.WTW.com

Xylem Analytics Germany GmbH
Dr.-Karl-Slevogt-Str. 1
82362 Weilheim
Germany