


# Prisma

## ETUVE MINI EM 4 Sans portoir



## MODE D'EMPLOI ETUVE MINI EM 4 Sans portoir

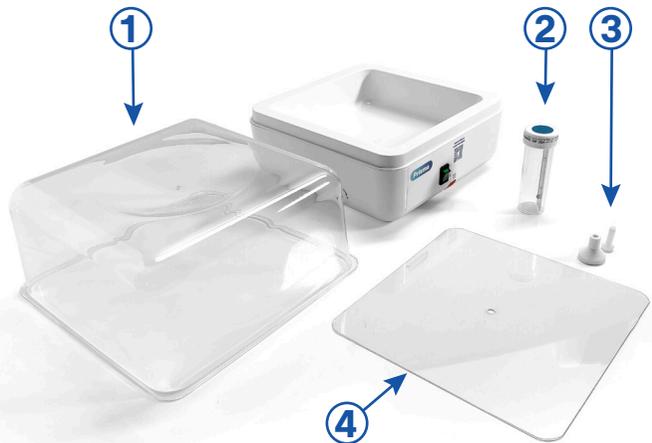

### COMPOSITION :

#### Étuve livrée avec :

- Une poignée (3) assemblant le couvercle (1) et le plateau (4)
- Un thermomètre de contrôle (2)

### CARACTÉRISTIQUES :

- Température idéale de la pièce : 20° C
- Réglage de la température de l'étuve (5) : de l'ambiance +5° C, jusqu'à 44° C
- Volume utile : 4 litres
- Dimensions externes : 23 x 23 x h. 18 cm
- Dimensions internes : 18 x 18 x h. 12,5 cm
- Poids : 1,5 kg
- Puissance : 25 W / 220 V.

### MISE EN MARCHÉ :

- Mettre sous tension avec l'interrupteur (6)
- Régler la température à l'aide de la vis (5)
- Lorsque l'étuve chauffe, le voyant lumineux rouge s'allume (7)
- Ne pas mettre d'eau dans l'étuve (sinon : exclusion de la garantie)

## DIRECTIONS FOR USE INCUBATOR MINI EM 4 Without holder



### COMPOSITION :

#### Incubator delivered with :

- One screwing button (2) assembling the lid (1) and the shelf (4)
- One control thermometer (2)

### FEATURES :

- Ideal temperature of the room : 20° C
- Operating range : from +5°C to 44°C with the help of the adjustment crew (5)
- Volume : 4 L.
- Overall dimensions : 23 x 23 x h. 18 cm
- Inner dimensions : 18 x 18 x h. 12,5 cm
- Weight : 1,5 kg
- Consumption : 25W / 220V

### FUNCTIONING :

- Switch on (6)
- Adjust the temperature with screw (5)
- Light (7) shows when incubator warms up
- Do not put water in the incubator (otherwise: warranty void)