LANGE 

DOC022.98.90414

LT20

12/2014, Edition 2

Основно ръководство за потребителя	bg
Grundlæggende brugervejledning	da
Basis Bedienungsanleitung	de
Βασικό Εγχειρίδιο Χρήστη	el
Basic User Manual	en
Manual básico del usuario	es
Basic käyttöohjeet	fi
Manuel d'utilisation de base	fr
Osnovni korisnički priručnik	hr
Alap felhasználói kézikönyv	hu
Manuale utente di base	it
Basisgebruikershandleiding	nl
Podstawowa instrukcja obsługi	pl
Manual de bază al utilizatorului	ro
Основное руководство пользователя	ru
Základný návod na použitie	sk
Osnovni uporabniški priročnik	sl
Grundläggande bruksanvisning	sv
Temel Kullanıcı Kılavuzu	tr

Английски	5
Dansk	11
Deutsch	17
Ελληνικά	23
English	29
Español	35
Suomi	41
Français	47
Hrvatski	53
Magyar	59
Italiano	65
Nederlands	71
Polski	77
Română	83
Русский	89
Slovenščina	95
Slovensky	101

Svenska	107
Türkçe	113

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Technical data

Subject to changes.

Performance specifications	Dimensions and technical data
Principle	Temperature control unit for 27 sample cells
Dimensions (W x H x D)	200 x 185 x 230 mm mm
Weight	4 kg
Operating temperature	+10°C – +30 °C, maximum 80 % relative humidity (without condensation formation)
Storage temperature	–20 °C – +60 °C, maximum 80 % relative humidity (without condensation formation)
Power consumption	65 VA
Mains connection	Input: 230 V +/- 10 % / 50 Hz
Fuse	T 315 mA L; 250 V (x 2)

General information

Safety notes

Read through the entire manual carefully before you unpack the device, set up or put into operation. Respect all danger and warning notes. Non-observance could lead to serious injury of the operator or to damage to the device.

To make sure that the protection equipment of the device is not impaired, this device may be used or installed in no other way than that described in this manual.

Danger notes in this manual

DANGER

Indicates a potentially or imminently dangerous situation that, if it is not avoided, leads to death or to serious injuries.

WARNING

Indicates a potentially or imminently dangerous situation that, if it is not avoided, can lead to death or to serious injuries.

CAUTION

Indicates a possible dangerous situation that can have minor or moderate injuries as the result.

NOTICE

Indicates a situation that, if it is not avoided, can lead to damage to the device. Information that particularly should be emphasized.

Note: Information that supplements aspects from the main text.

Warning labels

Observe all marks and labels that are attached to the device. Non-observance can result in personal injury or damage to the device.

	This symbol is a warning triangle. Follow all safety notes that follow this symbol to prevent possible injuries. If this symbol is located on the device, it refers to information in the operating and/or safety notes of the user manual.
	This symbol can be attached to a housing or a barrier in the product and shows that electric shock risk and/or the risk of a death through electric shock exists.
	Electrical equipment marked with this symbol may as of August 12, 2005 Europe-wide no longer be disposed of in unsorted house or industrial waste. According to valid provisions (EU Directive 2002/96/EC), from this point consumers in the EU must return old electrical devices to the manufacturer for disposal. This is free for the consumer. Note: You obtain instructions on the correct disposal of all (marked and not marked) electrical products that were supplied or manufactured by Hach Lange at your relevant Hach Lange sales office.

WARNING

The device may not be used in dangerous environments.

The manufacturer and its suppliers accept neither an express or indirect guarantee for use with high-risk activities.

Follow in addition to the local valid guidelines the following safety notes.

Safety notes for the regulatory use of the device:

- Do not operate the device near highly flammable substances, such as fuels, easily combustible chemicals and explosives.
- Do not operate the device near combustible gases, vapors or dust.
- Do not subject the device to strong vibrations or impacts.
- Do not open the device.
- If you do not use the device in accordance with regulations, this leads to loss of the guarantee.

Chemical and biological safety

DANGER

Potential danger in the event of contact with chemical/biological substances. Working with chemical samples, standards and reagents is linked with dangers. Make yourself familiar with the necessary safety procedures and the correct handling of the chemicals before the work and read and follow all relevant safety data sheets.

With normal operation of this device the use of chemicals that pose a risk health or of biologically harmful samples can be required.

- Before handling these substances observe all danger notes and safety information printed on the containers of the original solutions and in the safety data sheet.
- Dispose of all consumed solutions in accordance with the national regulations and laws.
- Select the type of protective equipment according to the concentration and quantity of the dangerous substance at the respective work place.

Startup

Change of the fuse

DANGER

Some circuits in this device are protected by fuses against overvoltages. For constant protection against risk of fire, replace these fuses only with fuses of the same type and power rating.

DANGER

Defective fuses generally indicate a problem in the device. If fuse errors repeatedly occur, contact the service department for instructions on the return procedure for repair purposes. Under no circumstances attempt to repair the device independently.

Operating environment

Observe the following points so that the device functions faultlessly and thus has a long service life.

- Place the device securely on an even surface. Push no objects under the device.
- The ambient temperature for transport, storage and operation must be +10 to +30 °C (+50 – +86 °F).

NOTICE

Protect the device against extreme temperatures from heaters, direct sunlight and other heat sources.

- The relative humidity must be under 80 % (no condensate formation).
- Keep above and on all sides of the device a clearance of at least 15 cm, so that air can circulate and overheating of the electrical parts is avoided.

- Do not use or store the device in extremely dusty, humid or wet places.

⚠ WARNING

Electrical dangers and fire hazard. Use only the supplied power cable.
Only qualified experts may perform the tasks described in this section of the manual, in compliance with all locally applicable safety regulations.

NOTICE


Use only an earthed socket for the connection of this device to the power supply.
If you are not sure if the socket is earthed, have this checked by a qualified electrician.

The power plug serves in addition to the power supply to isolate the device quickly from the mains where necessary.

This is recommended for long-term disuse and can prevent potential dangers in the event of a fault.

Therefore make sure that the socket to which the device is connected is easy to reach by each user at all times.

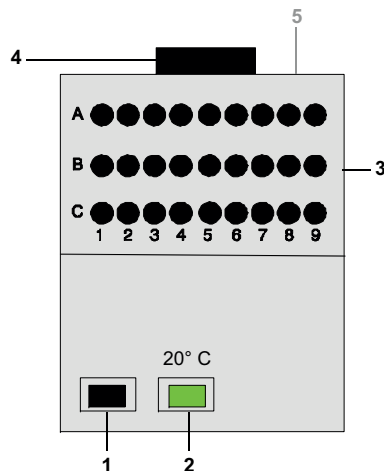Figure 1 LT20 equipment configuration

1	On/off switch	4	Fan
2	Temperature indicator	5	Rear: mains connection with fuse
3	Temperature control block (openings for 27 sample cells)		

The power cable is connected on the rear of the thermostat to an earthed mains socket (230 volt +/- 10 % / 50 Hz). The device is switched on with the toggle switch (on/off switch).

Note: Do not turn the device off and on in rapid succession. Always wait approximately 20 seconds before switching on again so as not to damage the electronics and mechanics of the device.

Routine operation

- Prepare the cells in accordance with the current operation procedure.
- Switch on the thermostat.
- Insert the cells in the temperature control block.
- The thermostat should only be operated when the cover is in place. If the green lamp goes out, the temperature is higher or lower than 20°C.

Clean the device

CAUTION

Potential dangers through pinches, burns and chemicals and for the eyes. Before all cleaning work, always disconnect the device from the power supply.

NOTICE

Under no circumstances may solvents like turpentine, acetone or similar be used to clean the device, including accessories.

Device

- Clean the housing and all accessories only with a soft, damp cloth. Mild soap solution can also be used.
- Allow no water to get into the temperature control block.
- Use cotton buds to clean or dry the cell housing.

Warranty and liability

The manufacturer warrants that the supplied product is free of material and manufacturing defects, and undertakes to repair or to replace any defective parts without charge.

The guarantee period for devices is 24 months. Upon conclusion of a maintenance contract within the first 6 months after purchase, the guarantee period extends to 60 months.

For defects to which the lack of assured properties also counts, the supplier is liable with exclusion of further claims as follows: All those parts that, within the guarantee period calculated from the day of the risk transfer, can be demonstrated as a result of circumstances prior to the risk transfer, in particular due to defective design, inferior materials or defective execution, to be unusable or whose usability is significantly impaired, are improved or replaced in accordance with the choice of the supplier free of charge. The identification of such defects must be notified to the supplier immediately, however at the latest seven days after the identification of the fault, in writing. If the customer omits this notification, the performance is considered approved despite the defect. Further liability for indirect or direct damages does not exist. If device-specific maintenance or inspection work prescribed by the supplier is to be performed within the guarantee period by the customer (maintenance) or by the supplier (inspection) and these requirements are not met, claims for damages that result from non-observance of these requirements are void.

Further claims, in particular for consequential damages, cannot be made valid.

Wear parts and damages that arise through improper handling, incorrect installation or non-regulatory use are excluded from this clause.

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ANALİZ SİSTEMLERİ LTD.ŞTİ.**