English

Español

Français

Deutsch

Italiano

Português

Svenska

Nederlands

Dansk

Polski

Čeština

Magyar

SF-110B Printer User Guide

Scan the QR code to download the full instruction manual or visit <http://www.ohaus.com/support> to view the full user manual.

1 INTRODUCTION

Thank you for purchasing this product.

This User Guide contains installation, operation, and maintenance instructions for SF110B printer. Please read it completely before installation and operation. For further detail information, consult the Instruction Manual on www.ohaus.com.

1.1 Safety Precautions

Signal Words

WARNING For a hazardous situation with medium risk, possibly resulting in severe injuries or death if not avoided.

CAUTION For a hazardous situation with low risk, resulting in damage to the device or the property or in loss of data, or minor or medium injuries

ATTENTION For important information about the product. May lead to equipment damage if not avoided.

NOTE For useful information about the product.

Warning Symbols

General hazard

Explosion hazard

Safety Precautions

CAUTION: Read all safety warnings before installing, making connections, or servicing this equipment. Failure to comply with these warnings could result in personal injury and/or property damage. Retain all instructions for future reference.

- Before connecting power, verify that the AC adapter's input voltage range and plug type are compatible with the local AC mains power supply.
- Do not position the equipment such that it is difficult to reach the power connection.
- Make sure that the power cord does not pose a potential obstacle or tripping hazard.
- Operate the equipment only under ambient conditions specified in these instructions.
- This equipment is for indoor use only.
- Do not operate the equipment in wet, hazardous, or unstable environments.
- Do not allow liquids to enter the equipment.
- Do not place the equipment upside down on the platform.
- Use only approved accessories and peripherals.
- Disconnect the equipment from the power supply when cleaning.
- Service should only be performed by authorized personnel.

WARNING: Never work in an environment subject to explosion hazards! The housing of the instrument is not gas tight. (Explosion hazard due to spark formation, corrosion caused by the ingress of gases)

1.2 Intended Use

This instrument is intended for use in businesses and light industry. It shall only be used for measuring the parameters described in these operating instructions. Any other type of use and operation beyond the limits of technical specifications, without written consent from OHAUS, is considered as not intended. This instrument complies with current industry standards and the recognized safety regulations; however, it can constitute a hazard in use.

If the instrument is not used according to these operating instructions, the intended protection provided by the instrument may be impaired.

2 INSTALLATION

2.1 Unpacking

Carefully remove your SF-110B printer and each of its components from the package. Save the packaging to ensure safe storage and transport.

Please read the manual completely before installing and using the printer to avoid incorrect operation. Components included:

- Printer
- User Guide
- Power Adapter with plugs*
- Paper Roll
- RS232 Cable
- Paper Rotary Axis
- Ink Ribbon

Note: * The included plugs vary depending on the region.

2.2 Connecting and Printing

Before connecting, make sure the instrument and printer have the same RS232 parameter setting. For instrument Baud rate settings, please see the related manual.

- Choose and install the proper power adapter converter according to the located country/region.
- Disconnect the power supply.
- Connect the printer to the instrument with the RS232 cable.
- Connect the printer to the mains power and use the main switch <I/O> to power on the printer.
- Press **PRINT** key on the printer or on the instrument.


Note: Depending on the output setting of the weighing instrument, it may be necessary to wait for the data to stabilize after pressing the print key, but not immediately.

3 OPERATION


3.1 Overview

3.1.1 Parts and Features

1. Paper cover
2. Paper Cutter
3. Control panel
4. Power Switch
5. Power jack
6. USB device interface
7. RS232 interface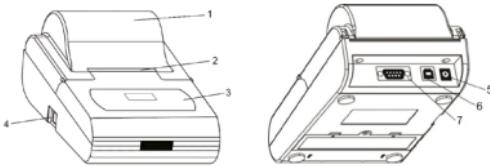

3.1.2 Control Panel

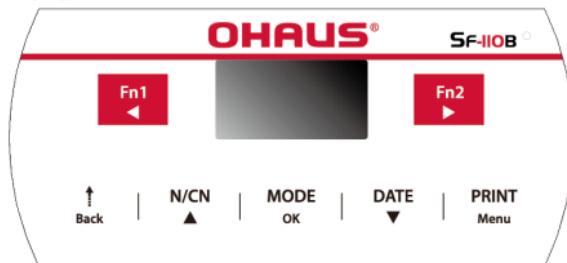Buttons	Functions
	Fn1 <ul style="list-style-type: none"> - Short press: Enter the sample name and lot number. - Press (in settings menu): Move the input position to the left.
	Fn2 <ul style="list-style-type: none"> - Short press: Open the related function menu of the current operation mode. - Long press: Enable or disable the automatic tare function. - Press (in settings menu): Move the input position to the right.
	Back <ul style="list-style-type: none"> - Short press: Feed the printer by one line. - Long press: Continuously feed until released. - Short press (in settings menu): Back to previous menu or cancel current parameter modification.

Buttons	Functions
N/CN ▲	N/CN <ul style="list-style-type: none"> - Short press: Enable or disable the numbering function. - Long press: Reset the number to "000". - Lone press (more than 4 seconds): modify the report's serial number. ▲ <ul style="list-style-type: none"> - Short press (in settings menu): Switch to the previous setting options or increase parameter value. - Long press (in settings menu): Quickly increase the value.
MODE OK	MODE <ul style="list-style-type: none"> - Short press: Choose operation modes. - Long press: Execute the user login / logout operation. OK <ul style="list-style-type: none"> - Short press (in settings menu): Access to a setting item or confirm the modified parameters.
DATE ▼	DATE <ul style="list-style-type: none"> - Short press: Print the current date and time. - Long press: Enter the Date/Time Settings menu. ▼ <ul style="list-style-type: none"> - Short press (in settings menu): Switch to the next setting item or decrease the parameters setting value. - Lone press (in settings menu): Quickly decrease the value.
PRINT Menu	PRINT <ul style="list-style-type: none"> - Short press: Perform printing operation. Menu <ul style="list-style-type: none"> - Long press: Enter the settings menu.

Notes:

Short Press: Press less than 2 seconds.

Long Press: Press and hold for more than 2 seconds.

3.1.3 Display

Status _____

Mode _____

Time _____

Status

The content of the status bar varies depending on the operation mode.

NAME	Sample name / lot number printout.
O/T	Enable Auto-tare.

Mode: Indicates the currently selected mode of operation.**Time:** Displays the current date and time. The current time is always displayed in most modes.

4 MAINTENANCE

Under normal conditions the printer requires practically no maintenance. The OHAUS service department is available if servicing is required. Please ask your responsible OHAUS dealer for details.

- When cleaning the printer, make sure no liquid flows into the printer.
- Keep the irritant liquid far away from printer.
- Remove the power cables and the data cables before cleaning.
- Please use soft dry cloth to wipe the printer after cleaning.
- Dry your hands before coming into contact with the AC adaptor.
- Keep it in a dry environment, if the printer will not be used for a long time pack it into the factory packaging.
- To avoid damage to the printer, it is strongly recommended to use ink ribbon and other consumables from the original supplier.
- Turn off the power supply of printer after usage. Please remove the adaptor/batteries if it is not going to be used for a long time.
- Always use original consumables. Unqualified consumables will damage the printer or shorten the life of the printer.

5 TECHNICAL DATA

5.1 Equipment Ratings

Indoor use only

Altitude:	2000m
Operating temperature:	0 to 40°C
Humidity:	Maximum relative humidity 80% for temperatures up to 31 °C decreasing linearly to 50% relative humidity at 40°C.
Electrical supply:	12VDC, 1A. (For use with certified or approved power supply, which must have a SELV and limited energy output.)
Voltage fluctuations:	Mains supply voltage fluctuations up to $\pm 10\%$ of the nominal voltage.
Oversupply category (Installation category):	II
Pollution degree:	2

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P/N 30719423 A © 2022 OHAUS Corporation, all rights reserved/ todos los derechos reservados / tous droits réservés / alle Rechte vorbehalten / tutti i diritti riservati / todos os direitos reservados / Alla rättigheter förbehållna / alle rechten voorbehouden / alle rettigheder forbeholdes / wszelkie prawa zastrzeżone / všechna práva vyhrazena / Minden jog fenntartva
OHAUS Corporation