Kit Gram-Hücker

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Kit Gram-Hücker

STAINING KIT FOR THE DIFFERENTIATION OF GRAM-POSITIVE FROM GRAM-NEGATIVE BACTERIA

REF Fast action : 361520-0000
361525-0000
Slow action : 362860-0000

Principe :

Gram-Hücker staining is a differential staining based on the permeability of the bacterial wall. In this technique, the bacterial wall is not stained but its structure permits classification of Gram-positive or Gram-negative bacteria. Lugol solution permits the formation of an intracellular complex with Crystal violet oxalate. A more important permeability of Gram-negative bacteria wall allows alcohol to eliminate this complex. Gram-negative bacteria can fix Safranin and then appear stained orange-pink. Gram-positive bacteria, characterized by a less important permeability of wall, are not discoloured by Alcohol and remain stained violet.

Kit description :

Bottle 1 : Cristal violet oxalate	1 x 240ml or 1 x 1000 ml
Bottle 2 : Lugol, PVP-stabilized solution	1 x 240ml or 1 x 1000 ml
Bottle 3 : Fast differentiator (alcohol/acetone)	1 x 240ml or 1 x 1000 ml
or	
Bottle 5 : Slow differentiator (alcohol-based)	1 x 240ml
Bottle 4 : Safranin	1 x 240ml or 1 x 1000 ml

Number of tests possible: 230 to 300 slides or 950 to 1250 slides.

Use:

For professional use only.

In vitro use only. **IVD**
The collection and processing of chemical and biological waste must be conducted by specialized and registered companies.
Storage temperature : 15-25°C away from light.
Opened bottle shelf life: report to expiration date on the label.

Specimen preparation :

Specimen must be treated in accordance with procedures available in the laboratory and promulgated by national authorities. It is necessary to perform a preliminary fixation by heat and/or hot alcohol (cf. Notice 03: Smears fixation, prior to staining in Bacteriology).

Recommendations and/or notes of use:

Adding of Polyvinylpyrrolidone (PVP) to Lugol solution helps to avoid iodine migration and then provides a satisfactory stability of the ready-to-use plastic bottle packaged solutions. One can improve the sharpness of the observation with a green light-microscope (a green filter or a yellow one superposed on the blue filter). The contrast depends on the way one carry out the differentiation. The genus *Campylobacter* is badly coloured by Safranin, while the genus *Legionella* remains colourless.

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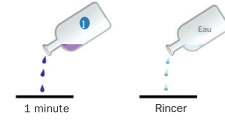
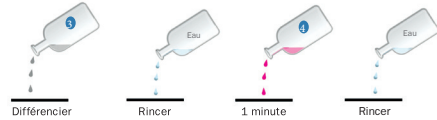
Kit Gram-Hücker

STAINING KIT FOR THE DIFFERENTIATION OF GRAM-POSITIVE FROM GRAM-NEGATIVE BACTERIA

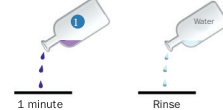
Staining procedure: Processing time : 3 minutes

Staining procedure:

Please read all this information carefully before using this device.

Processing time : 3 minutes

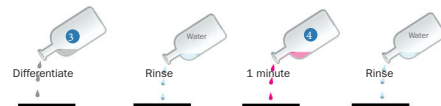
STEP 1 :
 • Place the slide on a stand with fixed smear on top.
 • Cover the smear with Crystal violet oxalate (bottle 1) for 1 minute.
 • Get rid of the stain and rinse with tap water to remove excess of Crystal violet oxalate.

STEP 2 :
 • Rinse with a jet of Lugol, PVP-stabilized solution (bottle 2) to remove excess of rinsing water.
 • Cover the slide with Lugol, PVP-stabilized solution (bottle 2) for 30 seconds to 1 minute.
 • Thoroughly rinse with tap water.

STEP 3 and 4 :
 • Discolour with the Fast differentiator (alcohol/acetone) for 2 to 5 sec or the Slow differentiator (alcohol-based) (bottle 5) for 20 to 40 sec.
 • Quickly rinse with tap water.
 • Cover the slide with Safranin (bottle 4) for 1 minute.
 • Quickly rinse with tap water.
 • Air dry.

Microscopic examination with a x100 immersion-objective.

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EN: DANGER. H225 Highly flammable liquid and vapour. H319 Causes serious eye irritation. H336 May cause drowsiness or dizziness. P210 Keep away from heat, hot surfaces, sparks, open flames and other ignition sources. No smoking. P261 Avoid breathing dust/fume/gas/mist/vapours/spray. P280 Wear protective gloves, protective clothing, eye protection. P312 Call a POISON CENTER or doctor/physician if you feel unwell. P337+P313 Eye irritation persists: Get medical advice/attention.

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EN: DANGER. H225 Highly flammable liquid and vapour. H318 Causes serious eye damage. P280 Wear protective gloves, protective clothing, eye protection. P210 Keep away from heat, hot surfaces, sparks, open flames and other ignition sources. No smoking. P305+P351+P338+P315 IF IN EYES: Rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing. Immediately call a POISON CENTER or doctor/physician.

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