Quanti-Tray*

Insert and Most Probable Number (MPN) Table

IDX 33/01 – 11/09
IDX 33/03 – 10/13
WATER ANALYSIS METHODS
www.afnor-validation.org

Quanti-Tray* Certificate of Sterility

This certifies that the enclosed Quanti-Trays have been sterilized with ethylene oxide.

For technical support, please call:

North/South America: +1 207 556 4496 or +1 800 321 0207

Europe: +00800 4339 9111

UK: +44 01638 676800

China: +86 21 61279528

Japan: +81 422 71 5921

Australia: +1300 44 33 99

idexx.com/water

Validated for use with IDEXX products only. Failure to use only IDEXX authorized products with this product will void any IDEXX warranty. IDEXX does not support any claims made by other manufacturers regarding product compatibility, performance or outcome.

The method Colilert*-18/Quanti-Tray* for water analysis is granted NF Validation by AFNOR Certification as an alternative method to the standard ISO 9308-1 for enumeration of *Escherichia coli* β -glucuronidase positive and coliform bacteria β -galactosidase positive in drinking water (except bottled water), under the Certificate number: IDX 33/01 – 11/09. For more information about end of validity, please refer to the certificate NF Validation available on website mentioned above.

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The method Enterolert-DW*/Quanti-Tray* for water analysis is granted NF Validation by AFNOR Certification as an alternative method to the standard ISO 7899-2 for enumeration of intestinal enterococci in drinking water (except bottled water), under the Certificate number: IDX 33/03 – 10/13. For more information about end of validity, please refer to the certificate NF Validation available on website mentioned above.

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IDEXX

One IDEXX Drive
Westbrook, Maine 04092 USA

*Quanti-Tray and Colilert are trademarks or registered trademarks of IDEXX Laboratories, Inc. or its affiliates in the United States and/or other countries.

Patent information: idexx.com/patents.

© 2021 IDEXX Laboratories, Inc. All rights reserved. • 06-02030-18

Quanti-Tray*

English

Introduction




IDEXX Quanti-Tray* (98-14855-00, 98-21378-00) is designed to give quantitated bacterial counts of 100 mL samples using IDEXX reagent products. Add the reagent/sample mixture to a Quanti-Tray, seal it in a Quanti-Tray* Sealer and incubate per the reagent instructions. Count the number of positive wells and use the Most Probable Number (MPN) Table attached to determine the MPN.

Contents



This package contains sterile 51-well Quanti-Trays.

User Instructions

1. Use one hand to hold a Quanti-Tray upright with the well side facing the palm.
2. Squeeze the upper part of the Quanti-Tray so that the Quanti-Tray bends toward the palm.
3. Open the Quanti-Tray by pulling the foil tab away from the well side. Avoid touching the inside of the foil or tray.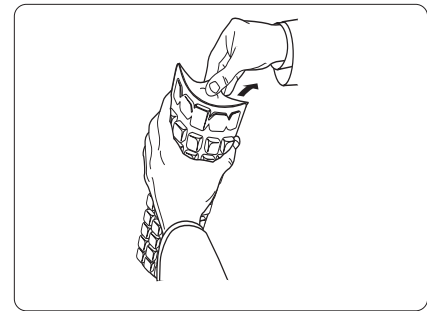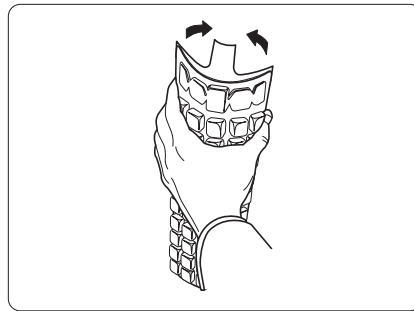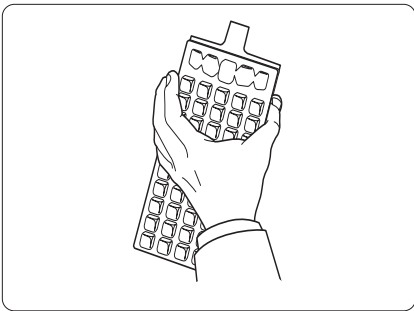

4. Pour the reagent/sample mixture directly into the Quanti-Tray, avoiding contact with the foil tab. Allow foam to settle.
5. Place the sample-filled Quanti-Tray onto the rubber tray carrier of the Quanti-Tray Sealer with the well side (plastic) of the Quanti-Tray facing down to fit into the carrier.
6. Seal according to the Quanti-Tray Sealer instructions.
7. Incubate according to reagent instructions.

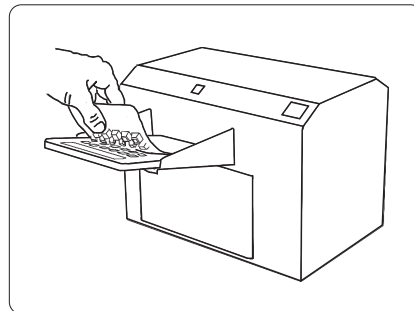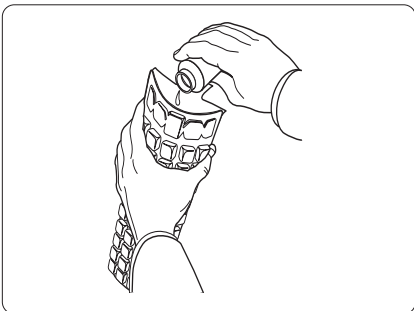8. Count positive wells and refer to the MPN table on the back of this instruction sheet to find the MPN.[†]
9. Dispose of media in accordance with good laboratory practices.

For technical support, please call:

North/South America: +1 207 556 4496 or +1 800 321 0207

Europe: +00800 4339 9111

UK: +44 01638 676800

China: +86 21 61279528

Japan: +81 422 71 5921

Australia: +1300 44 33 99

idexx.com/water

[†]Download the IDEXX MPN generator software for automated Quanti-Tray results at idexx.com/mpngenerator.

*Quanti-Tray is a trademark or registered trademark of IDEXX Laboratories, Inc. in the United States and/or other countries.

© 2021 IDEXX Laboratories, Inc. All rights reserved.

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