SFX/STEREO Series

INSTRUCTION MANUAL

Model
MS-2
STX
SFX-31
SFX-32
SFX-33
SFX-34
SFX-51
SFX-52
SFX-91
SFX-91D
ST-30FX
ST-50Led

Ver. 2.0 2019

Table of Contents

1. Warning	3
2. Symbols and conventions	3
3. Safety Information	3
4. Intended use	3
5. Overview	4
5.1 MS-2	4
5.2 STX	5
5.3 SFX-31	6
5.4 SFX-32	7
5.5 SFX-33	8
5.6 SFX-34	9
5.7 SFX-51	10
5.8 SFX-52	11
5.9 SFX-91	12
5.10 SFX-91D	13
5.11 ST-30FX	14
5.12 ST-50Led	15
6. Unpacking	16
7. Using the microscope	16
7.1 Place the specimen on the stage	16
7.2 Turning on the light	16
7.3 Adjust interpupillary distance	16
7.4 Focus	17
7.5 Dioptic compensation	17
7.6 Magnification	17
7.7 Contrast disc	17
7.8 Inserting the batteries	18
7.9 Charge Battery LED	18
8. Installation of camera drivers and software	19
9. Initial settings	19
10. Micrometric Slide M-005	21
11. Maintenance	22
Equipment disposal	23

1. Warning

This microscope is a scientific precision instrument designed to last for many years with a minimum of maintenance. It is built to high optical and mechanical standards and to withstand daily use. We remind you that this manual contains important information on safety and maintenance, and that it must therefore be made accessible to the instrument users. We decline any responsibility deriving from incorrect instrument use that does not comply with this manual.

2. Symbols and conventions

The following chart is an illustrated glossary of the symbols that are used in this manual.

CAUTION

This symbol indicates a potential risk and alerts you to proceed with caution.

ELECTRICAL SHOCK

This symbol indicates a risk of electrical shock.

3. Safety Information

Avoiding Electrical Shock

Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off position. Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users have full responsibility to use this equipment safely. Please follow the guidelines below, and read this manual in its entirety to ensure safe operation of the unit.

4. Intended use


For teaching use only. Not intended for any animal or human therapeutic or diagnostic use.

5. Overview


5.1 MS-2


5.2 STX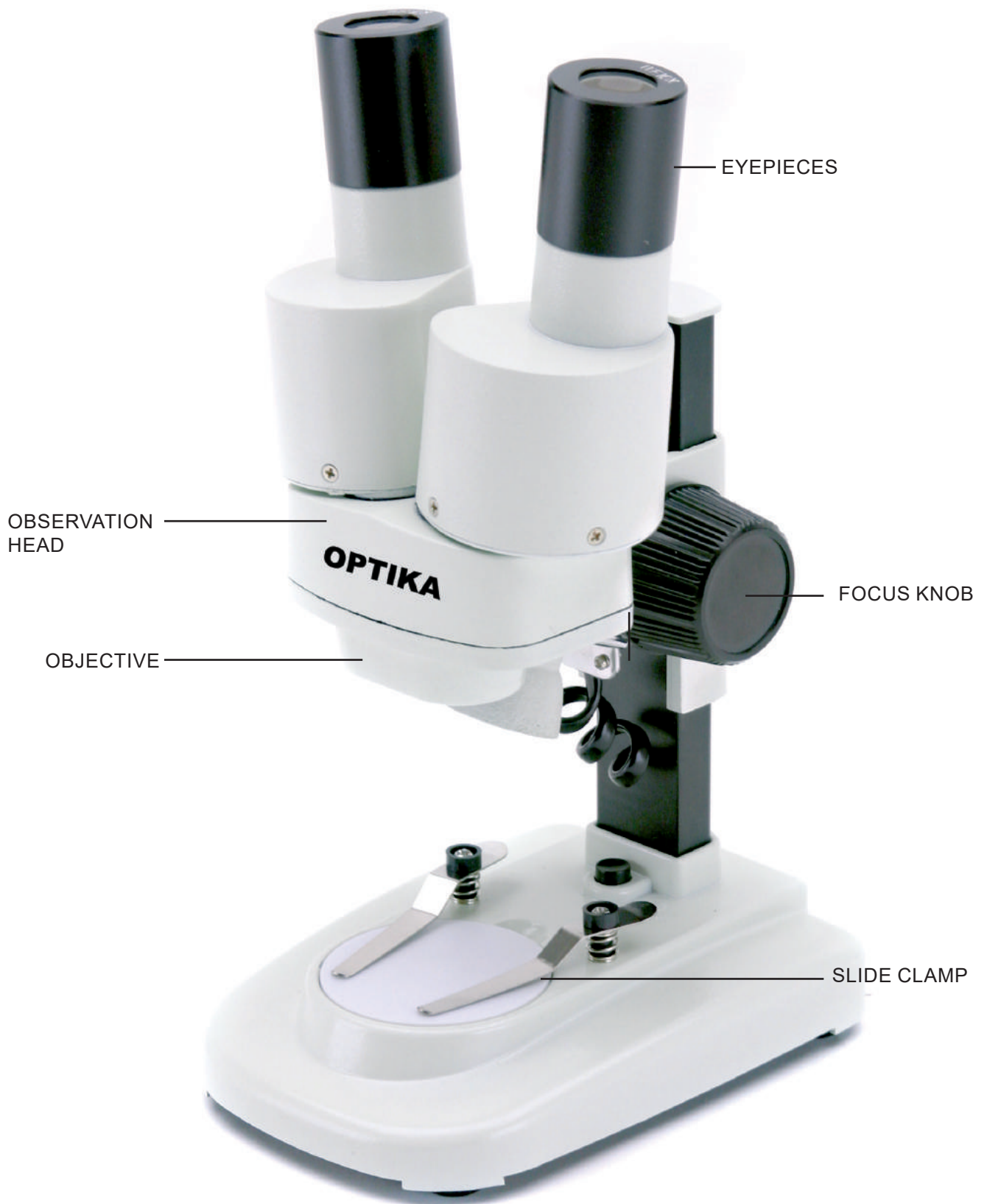


5.3 SFX-31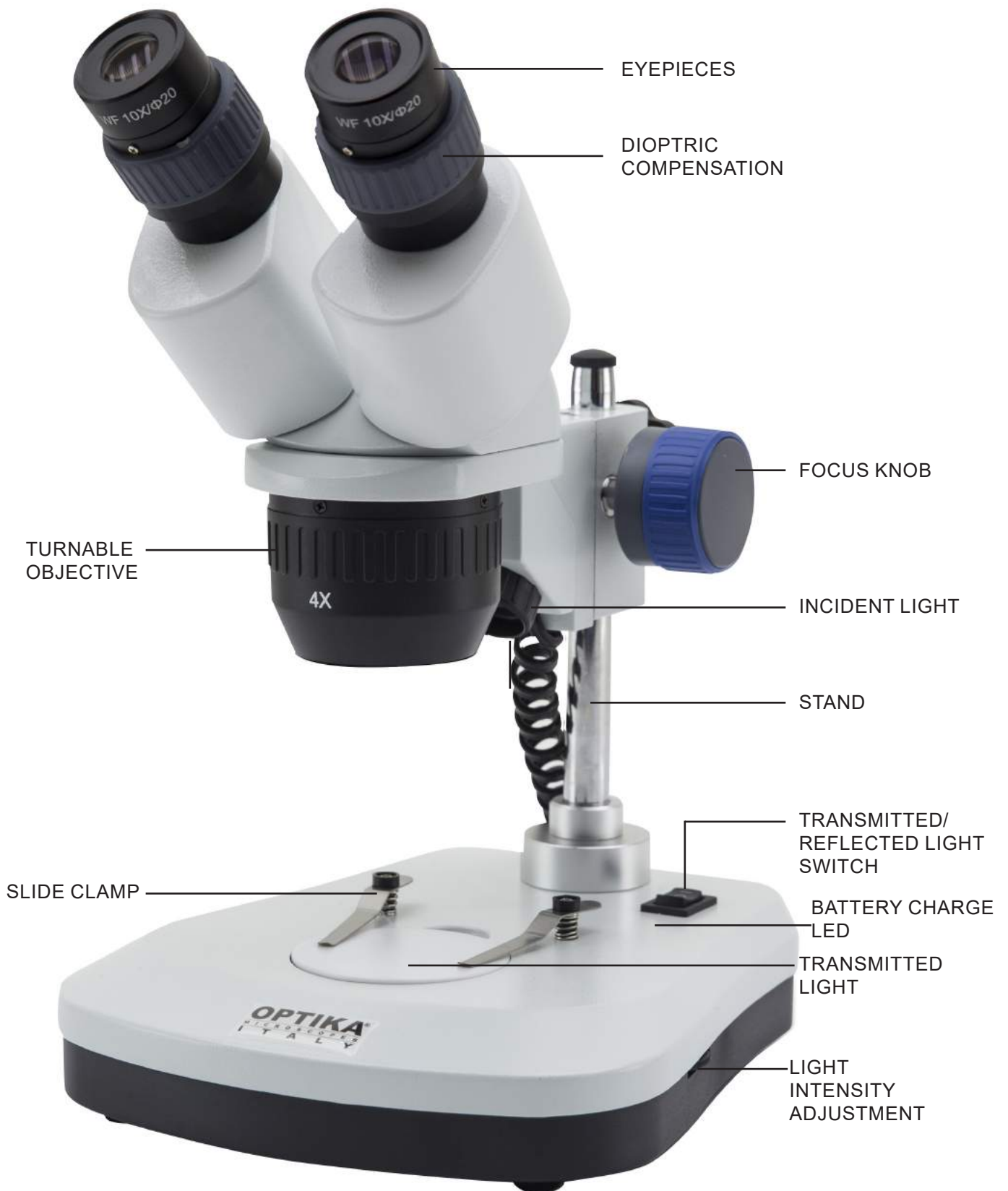


5.4 SFX-32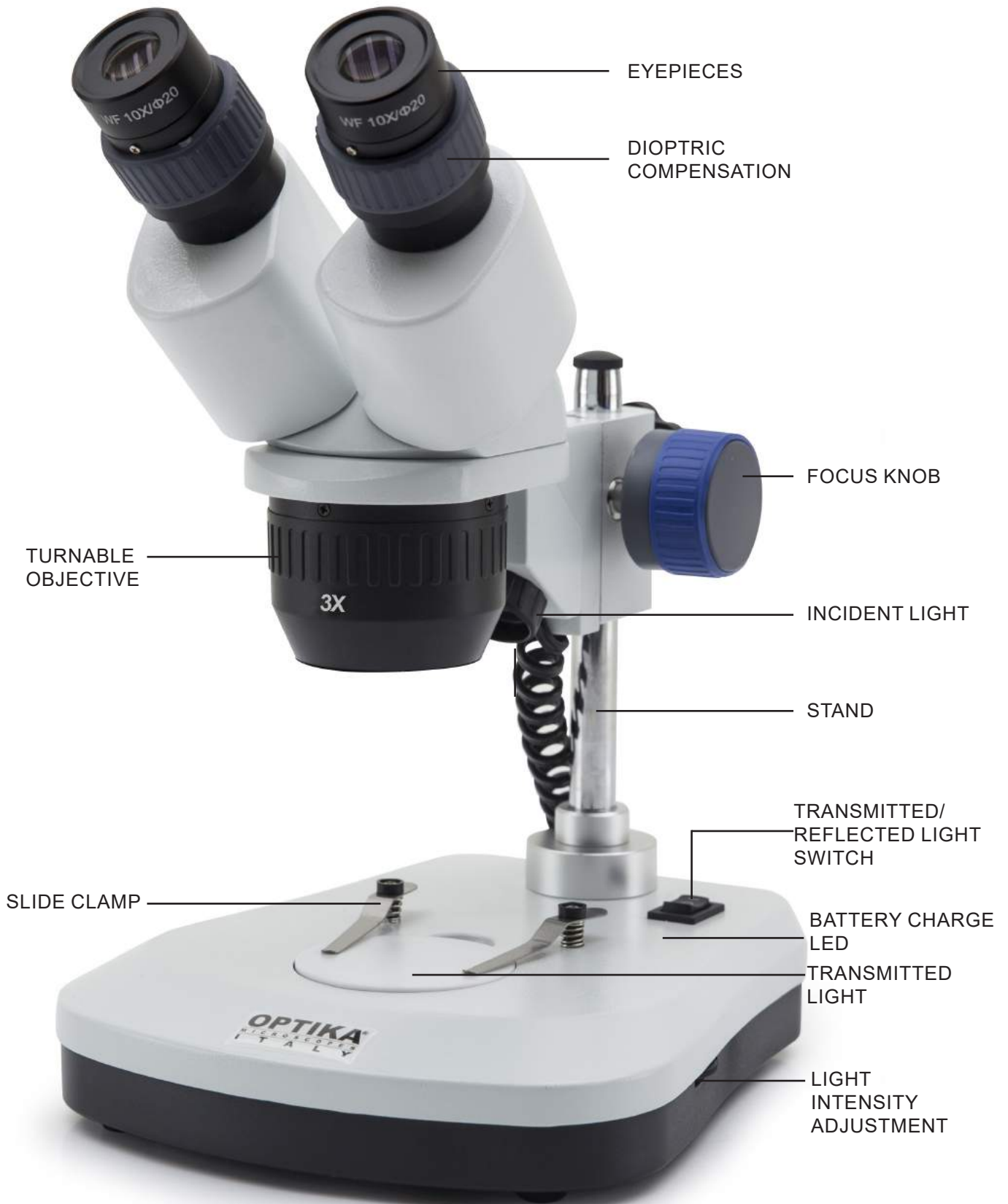

5.5 SFX-33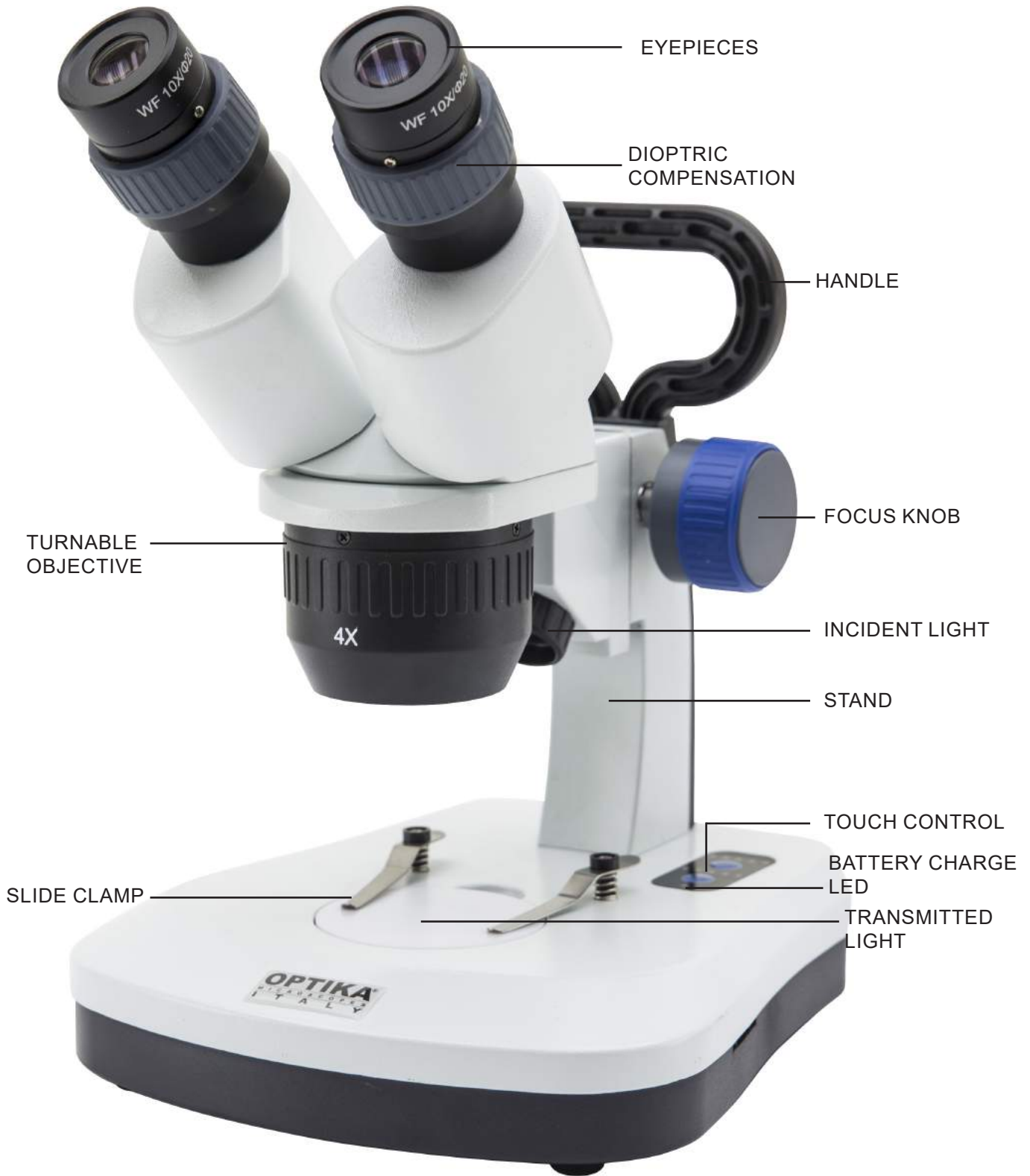

5.6 SFX-34

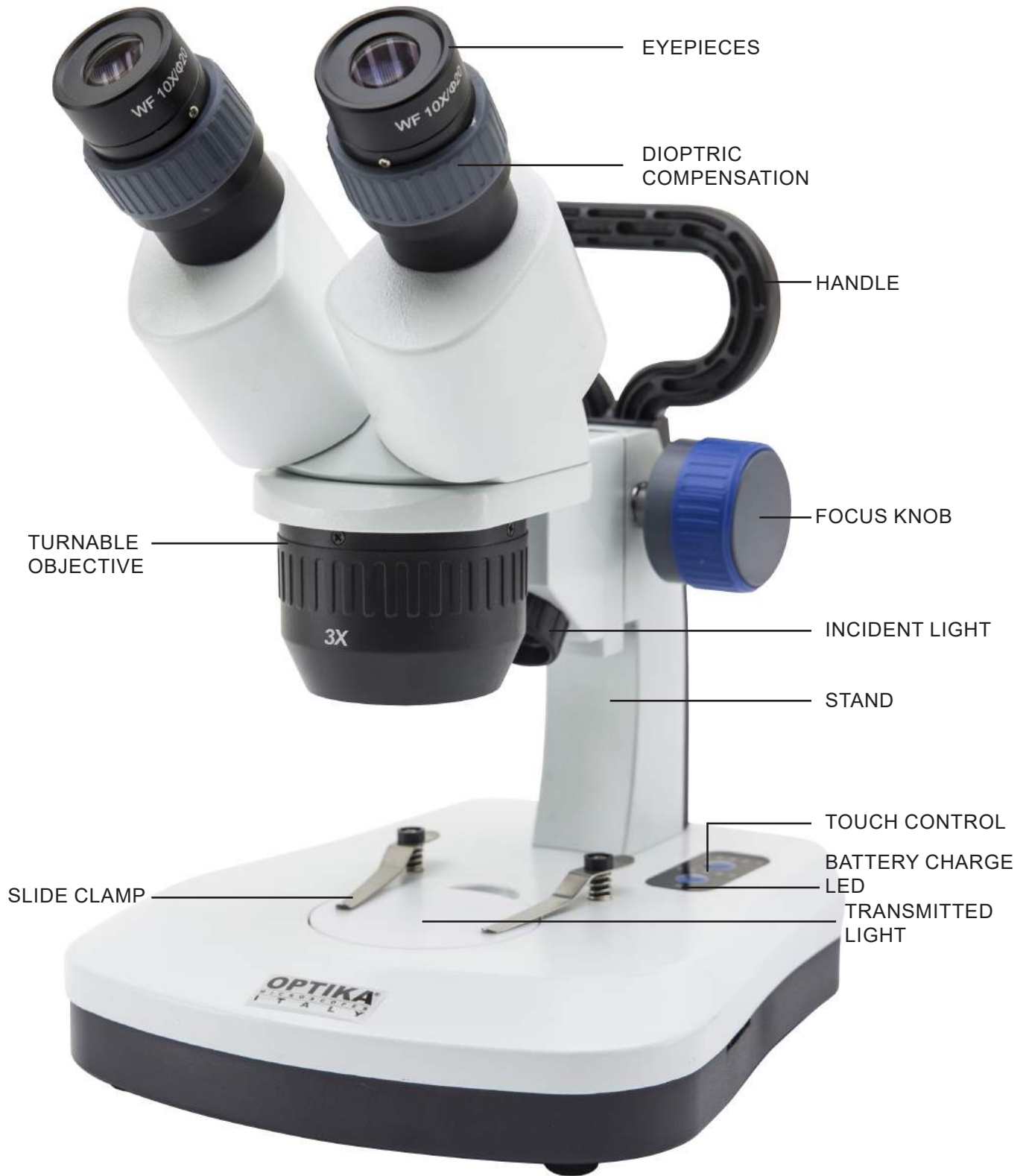
5.7 SFX-51


5.8 SFX-52


5.9 SFX-91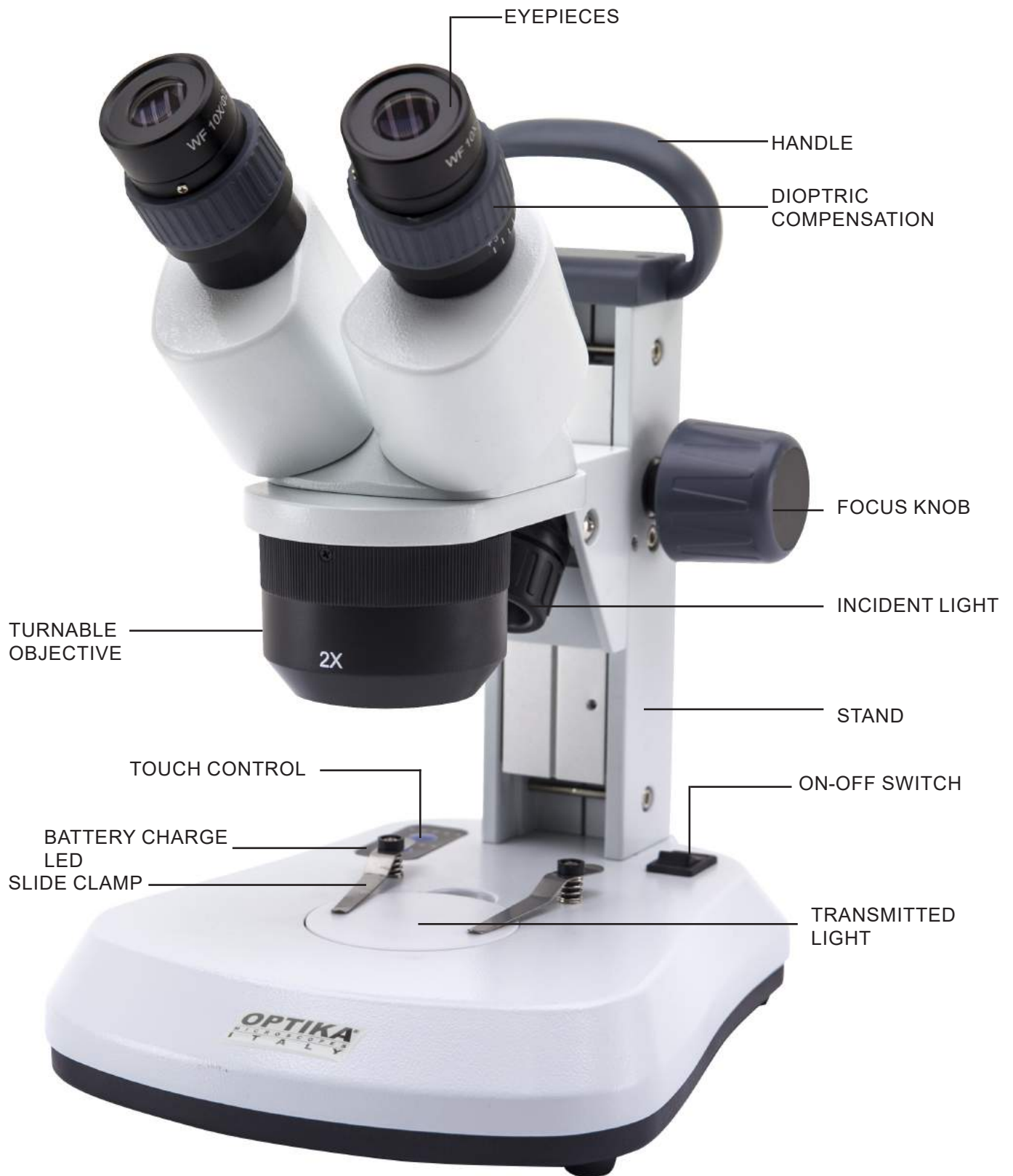


5.10 SFX-91D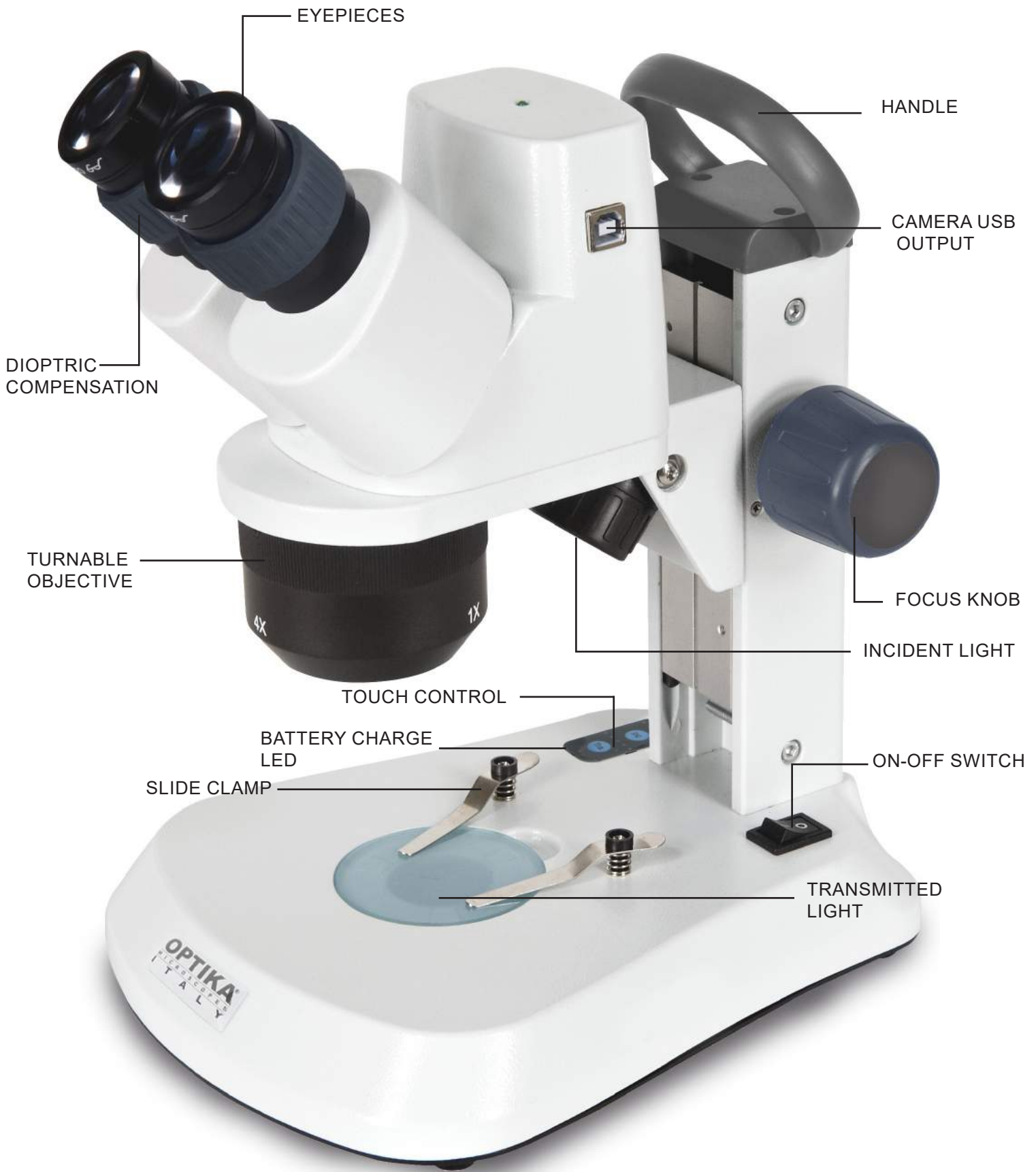

5.11 ST-30FX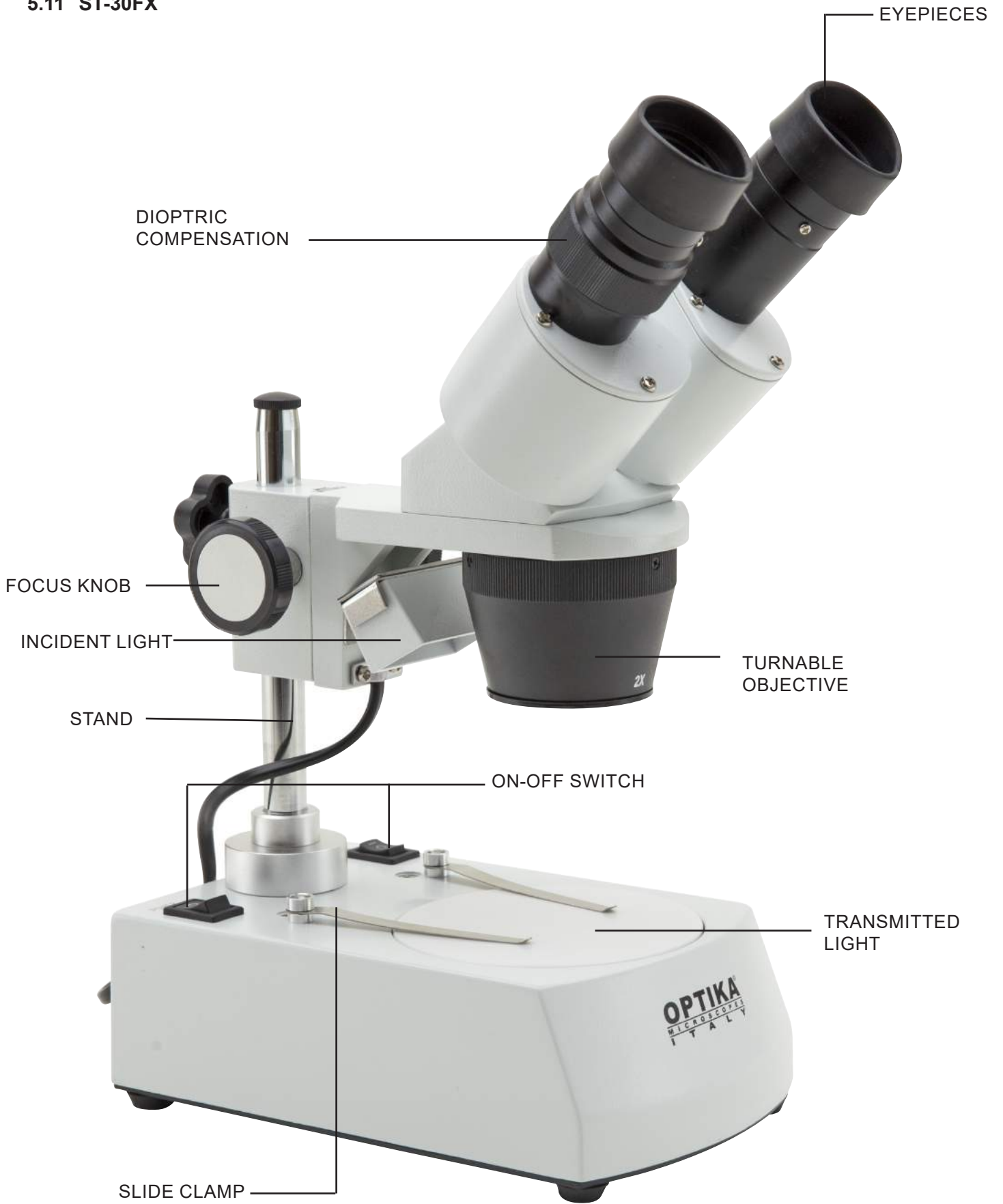

5.12 ST-50Led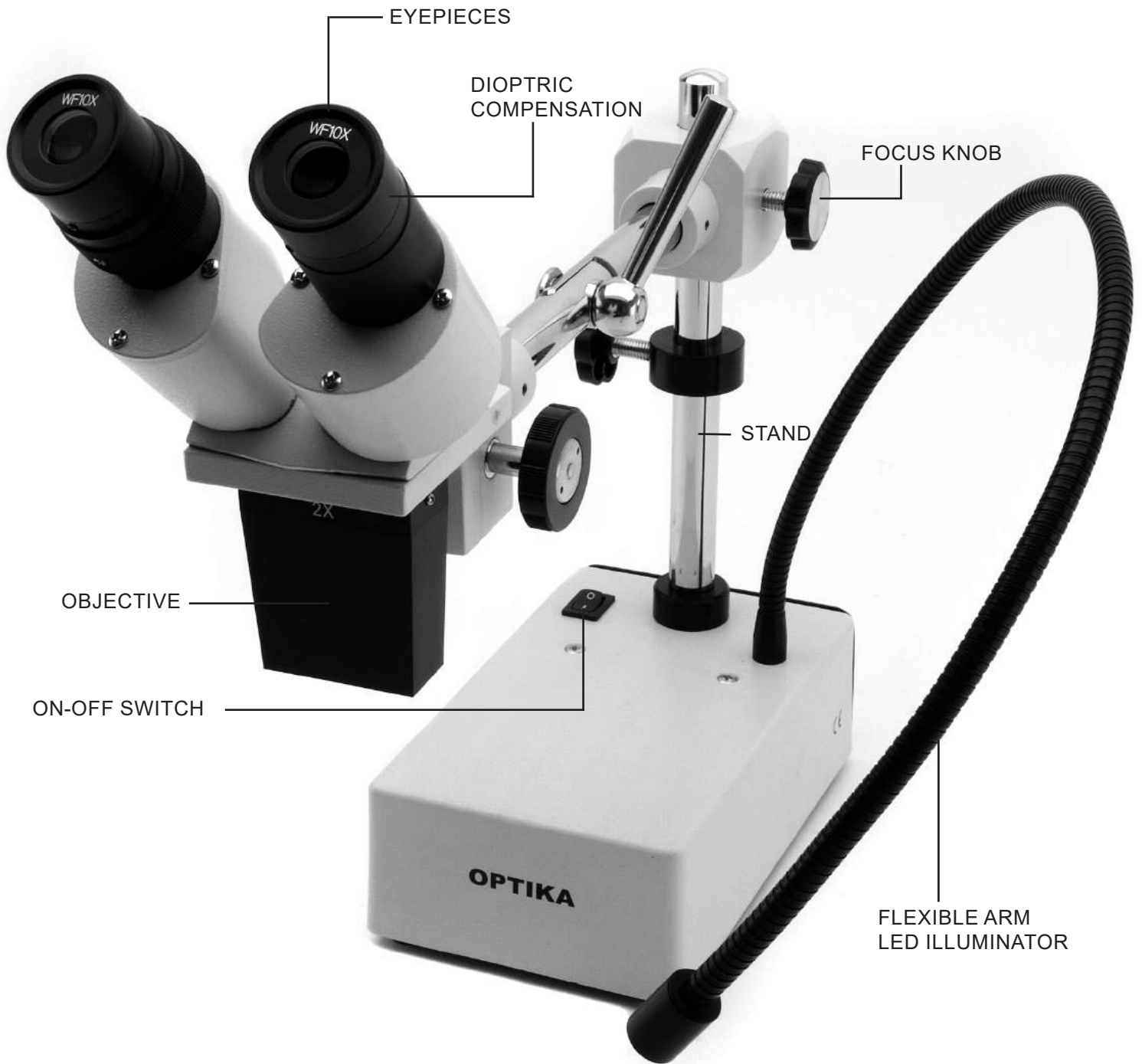

6. Unpacking

The microscope is housed in a moulded Styrofoam container. Remove the tape from the edge of the container and lift the top half of the container. Take some care to avoid that the optical items (objectives and eyepieces) fall out and get damaged. Using both hands (one around the arm and one around the base), lift the microscope from the container and put it on a stable desk.

7. Using the microscope

7.1 Place the specimen on the stage

Place the specimen on the microscope stage and lock the specimen using the slide-clamp if it is necessary. Ensure that the specimen is centred over the stage opening. (Fig.1)


7.2 Turning on the light

The microscope comes with an electrical illuminator. Insert the plug of the cable into the power socket (or the power supply jack in the socket), turn on the switch on the main body and select your light source. (Fig. 2A-2B)


For SFX models, repeatedly press the touch button in order to change the light intensity. Depending on the model, the light intensity can be adjusted with a knob on the right side of the stand. (Fig. 3-4)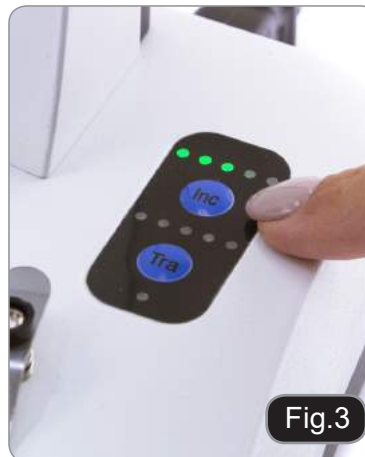

7.3 Adjust interpupillary distance

Hold the right and left parts of the observation head by both hands and adjust the interpupillary distance by turning the two parts until one circle of light can be seen. If two circles appear, the interpupillary distance is too big, and if two overlapped circles appear, the interpupillary distance is too small. (Fig. 5)

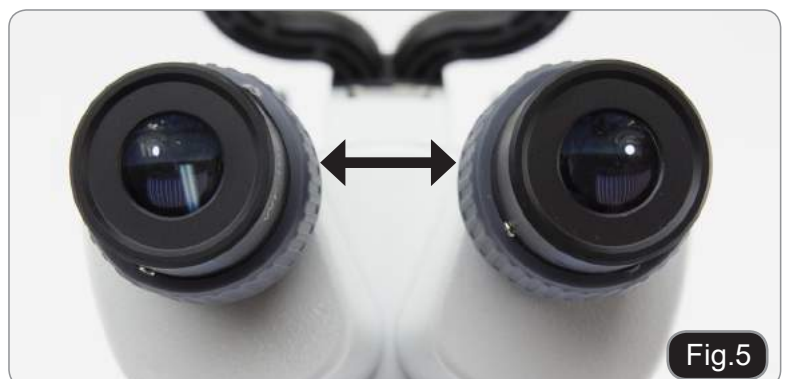7.4 Focus

Rotate the focusing knob to bring the sample into focus with the highest magnification. (Fig. 6)

Fig.6

7.5 Dioptic compensation

This compensation makes it possible for people with glasses to adjust the microscope to their eyes and use the microscope without glasses. Adjust the diopter compensation ring of the right eyepiece until the image of the right eyepiece is clear and sharp. Repeat the procedure for the left eyepiece. (Fig. 7)

Fig.7

7.6 Magnification

Select the desired magnification by rotating the objective. (Fig. 8)
Total magnification used can be calculated as:
Eyepiece magnification x Zoom magnification
x Objective lens magnification.

Fig.8

7.7 Contrast disc

You can use the black/white disc in order to enhance the contrast of the image when using the incident illumination. If you want to look at bright opaque objects, place the disc with the black side up. (Fig. 9)

Fig.9

7.8 Inserting the batteries

(Models with rechargeable batteries)

Before you start using the instrument, you must insert the rechargeable batteries.

Open the battery cover (located on the bottom of the instrument) and insert the provided batteries, taking care to respect the polarities. (Fig. 10-11)

When finished, close the battery cover again.

Fig.10

Fig.11

7.9 Charge Battery LED

(Only SFX models)

This LED ① shows the actual charge level of the batteries. (Fig. 12)

LED ON (green): batteries charged

LED OFF (no color): batteries to be recharged.

The status of the LED is independent of whether the power supply unit is connected or not.

Fig.12

8. Installation of camera drivers and software

Only for SFX-91D

Operating system requirements: Windows XP, Windows 7, Windows 8 or Windows 10.

If your PC is Windows XP: before the installation of your Camera please check if Service Pack2 is installed. If not, please update your XP from the Microsoft web site.


If your PC has a USB port less than 2.0, there may be some problems in capturing image from the camera due to the slow transmission.

1. Please insert your CD and run: `\\optika vision lite \ setup.exe`
2. Now Optika Vision Lite icon appears on your desktop.
3. Open the folder "**Drivers**" and run all the ".exe" files.
4. The installer may show warnings about missing Windows certification: ignore such warnings and select "**Continue**".
 - **Note : do NOT cancel the installation.**
5. At the end of the installation process restart the PC if requested.
6. Plug the Camera in your PC: the camera will be identified and Windows will recognize the driver which has just been installed from the CD.
Then ignore any message about Windows compatibility and select "**Continue**".
 - **Note : do NOT cancel the installation.**
7. System is ready.

9. Initial settings


Once started the software, the user must set some parameters in order to achieve a proper function of the whole system.

1. Click on the microscope icon. The following dialog appears: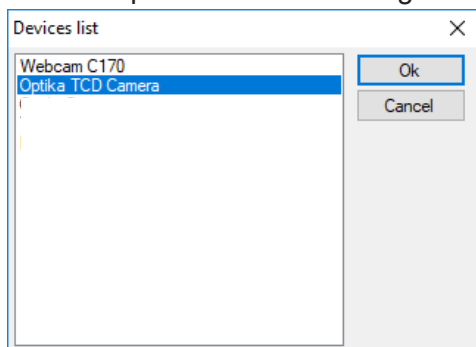



2. Select Optika TCD Camera.

3. A new dialog appears: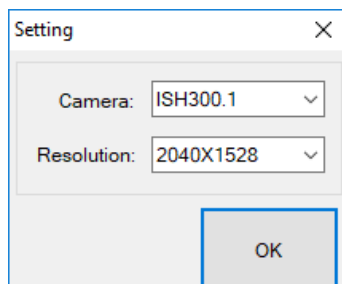

4. Press OK.

5. The following dialog showing the camera parameters appears: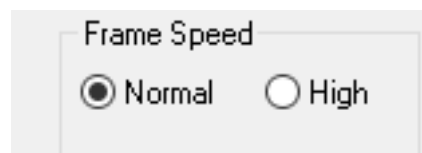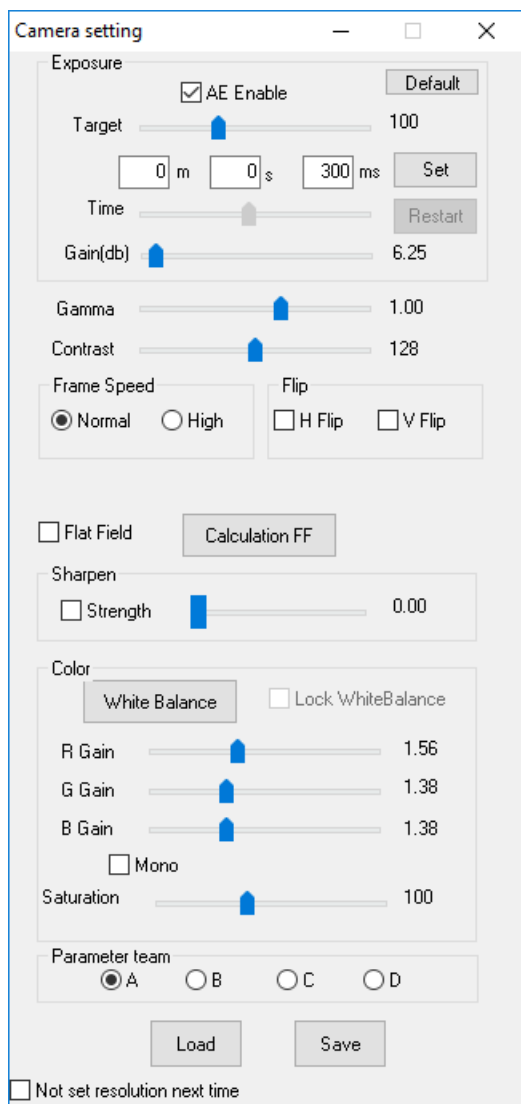


- **NOTE: if the user sets the Frame Speed to “High” all the software operations will dramatically drop down. We suggest to keep the setting to “Normal”.**

Please refer to the Optika Vision Lite instruction manual for the complete explanation of the software functions.


10. Micrometric Slide M-005

**Micrometric slide, 26x76mm, with 2 scales
(1mm/100div. for biological microscopes / 10mm/100div. for stereomicroscopes)**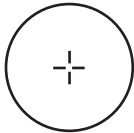

1 DIV=0.01mm

For biological microscopes calibration

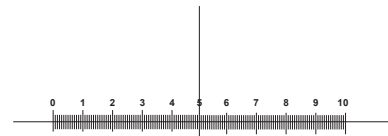1 DIV=0.1mm

For stereo microscopes calibration

11. Maintenance

Microscopy environment

This microscope is recommended to be used in a clean, dry and shock free environment with a temperature of 5°-40°C and a maximum relative humidity of 75 % (non condensing). Use a dehumidifier if needed.

To think about when and after using the microscope

- The microscope should always be kept vertically when moving it and be careful so that no moving parts, such as the eyepieces, fall out.
- Never mishandle or impose unnecessary force on the microscope.
- Never attempt to service the microscope yourself.
- After use, turn off the light immediately, cover the microscope with the included dust-cover, and keep it in a dry and clean place.

Electrical safety precautions

- Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off position.
- Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users do have full responsibility to use this equipment safely.

Cleaning the optics

- If the optical parts need to be cleaned try first to: use compressed air.
- If that is not sufficient: use a soft lint-free piece of cloth with water and a mild detergent.
- And as a final option: use the piece of cloth moistened with a 3:7 mixture of ethanol and ether.
- Note: ethanol and ether are highly flammable liquids. Do not use them near a heat source, near sparks or near electric equipment. Use these chemicals in a well ventilated room.
- Remember to never wipe the surface of any optical items with your hands. Fingerprints can damage the optics.
- Do not disassemble objectives or eyepieces in attempt to clean them.

For the best results, use the OPTIKA cleaning kit (see catalogue).

If you need to send the microscope to Optika for maintenance, please use the original packaging.

Equipment disposal

Art.13 Dlsg 25 July 2005 N°151. “According to directives 2002/95/EC, 2002/96/EC and 2003/108/EC relating to the reduction in the use of hazardous substances in electrical and electronic equipment and waste disposal.”

The basket symbol on equipment or on its box indicates that the product at the end of its useful life should be collected separately from other waste. The separate collection of this equipment at the end of its lifetime is organized and managed by the producer. The user will have to contact the manufacturer and follow the rules that he adopted for end-of-life equipment collection. The collection of the equipment for recycling, treatment and environmentally compatible disposal, helps to prevent possible adverse effects on the environment and health and promotes reuse and/or recycling of materials of the equipment. Improper disposal of the product involves the application of administrative penalties as provided by the laws in force.

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OPTIKA® S.r.l.