Bibby Sterilin



Assembly Instructions

- 1 Fasten the 4 hinge bars to panels using x2 M4 screws to each hinge component. Do not tighten at this stage.
- (2) Attach the handle to front door using x5 M4 screws.
- **3** Screw the countersunk metal screw and washer into the bottom of the front panel.
- 4 Connect the 2 door clips to the front of the side panels with x2 M3 screws and x2 dome nuts for each clip.
- (5) Fasten the tie bar with x2 M4 screws at each end. The unit should now be free standing with the tie bar supporting the 2 end units.
- (6) Insert the hinge bars connecting the front and back panels.
- 7 Squeeze some silicon compound into the grooves on the back and top panel so when assembled a good seal is formed.
- 8 Insert two side panels into back panel. Then place top panel into position.

List of components

x1 control panel (right x4 hinges hand side) x2 hinge bars x1 door panel x4 plugs x1 left panel x2 door clips

x1 top panel x2 door clips x1 top panel x1 magnet catcher

x1 back panel x1 countersunk screw & washer

x1 tie bar x4 dome nuts x1 door handle x4 M3 screws x1 tube of silicon rubber compound x17 M4 screws

