



LETHEEN SUPPLEMENT

Emulsione sterile di Lecitina per Letheen Agar Base e Letheen Broth Base
Sterile emulsion of Lecithin for Letheen Agar Base and Letheen Broth Base.

MODALITA' D'USO

Agitare bene il flacone e prelevare le seguenti quantità:
1. 10 ml da aggiungere ad 1 litro di Letheen Agar Base(codice 610076-620076); oppure
2. 7 ml da aggiungere ad 1 litro di Letheen Broth Base (codice 610208-620208)
Miscelare bene. Sterilizzare a 121 °C per 15 minuti.
Raffreddare a 50 gradi e distribuire in piastre petri o contenitori finali sterili.

INSTRUCTION FOR USE

Mix well the bottle and pick up the following quantities:
1. *10 ml to adding to 1 litre of Letheen Agar Base (Code 610076-620076); or*
2. *7 ml to adding to 1 litre of Letheen Broth Base (Code 610208-620208)*
Mix well. Sterilize at 121°C for 15 minutes.
Cool to 50°C and dispense in petri dishes or sterile final containers.

FORMULA (grammi per litro)

Lecitina..... 100 g
Acqua..... 1000 ml

FORMULA (grams per litre)

Lecithin..... 100 g
Water..... 1000 ml

CONTENUTO

4 Fiale x 50 ml di Letheen Supplement

CONTENT

4 Vials x 50 ml of Letheen Supplement

CONSERVAZIONE

2-8°C

STORE AT

5-12°C

VALIDITA'

2 anni

SHELF LIFE

2 years

PRODOTTO / PRODUCT	CODICE / CODE	CONFEZIONE / PACKAGING
LETHEEN SUPPLEMENT	80008	4 Fiale x 50 ml per *20 litri di Letheen Agar *28 litri di Letheen Broth 4 Vials x 50 ml for *20 liters of Letheen Agar *28 liters of Letheen Broth
LETHEEN AGAR BASE	610076	500g
LETHEEN AGAR BASE	620076	100 g
LETHEEN BROTH BASE	610208	500 g
LETHEEN BROTH BASE	620208	100 g

BIBLIOGRAPHY

1. Quisno, R., I. W. Gibby, and M.J.Foter. 1946. "A neutralizing medium for evaluating the germicidal potency of the quaternary ammonium salts. *Am.J.Pharm.* 118:320-323

Rev.01/01.02.2007