

Technical Data Sheet

Readycult®

Coliforms 50

Ordering number: 1.01295.0001

Selective enrichment broth for the simultaneous detection of total **coliforms** and **E. coli** within the bacteriological water examination.

General Information

Content: 20 snap packs

1 snap pack each for 50 ml of water sample.

Mode of Action

The high nutritional quality of the peptones and the incorporated phosphate buffer guarantee rapid growth of coliforms whereas lauryl sulfate largely inhibits the accompanying flora, especially the Gram-positive. By adding the chromogenic substrate X-GAL which is cleaved by coliforms and the fluorogenic substrate MUG which is highly specific for E. coli the simultaneous detection of total coliforms and E. coli is possible. The presence of total coliforms is indicated by a blue-green colour of the broth and E. coli by a blue fluorescence under UV-light.

Typical Composition (g/L)

Readycult® Coliforms 50	
Tryptose	0.25
Sodium chloride	0.25
Sorbitol	0.05
Tryptophan	0.05
di-Potassium hydrogen phosphate	0.135
Potassium dihydrogen phosphate	0.1
Lauryl sulfate sodium salt	0.005
X-GAL	0.004
MUG	0.0025
IPTG	0.005

pH: 6.8 ± 0.2 at 25°C

Preparation

1. Add 50 ml of water sample into a sterile, transparent vessel with screw cap. (minimum capacity: >50 ml) **Attention: please use material e.g. glass that is not self-fluorescing!**
2. Take one snap pack, shortly tap to ensure the granules are at the bottom. Bend the upper part of the snap pack until it breaks open. **Attention: do not touch the opening to avoid contamination risk!**
3. Add the content to the water sample. Seal the vessel and shake to dissolve the granules completely.

The prepared broth is clear and yellowish.

Incubation: up to 24 h at 35 °C to 37 °C.

If incubated at room temperature (20 to 25 °C) the incubation time is prolonged to 48 hours.

Experimental Procedure and Evaluation

Interpretation of results for the detection of Total Coliforms / E. coli:

Negative: No colour change.
The broth remains yellowish in colour.

Total coliforms: Any colour change of the broth to blue-green, even in the upper section of the broth only, confirms the presence of coliforms (X-GAL reaction).

No decolouration with shaking!

E. coli: Check blue-green coloured vessels for fluorescence by using UV-lamp (366 nm) in front of the vessel. Light blue fluorescence indicates presence of E. coli (MUG reaction).

Attention: Protect your eyes from direct UV light!

To confirm **E. coli** in the vessel with positive fluorescence, overlay the broth with 2.5 ml of KOVAC's reagent (indole reaction). A red ring confirms presence of **E. coli**.

	Colour change to blue-green	Fluorescence	Indole-Reaction
Total coliforms	+	-	-
E. coli	+	+	+
Negative	yellow colour (no change)		

Disposal

Autoclave the broth (15 min/121°C). Alternatively heat the broth for 30 min. in boiling water or use a proper disinfectant.

Storage

In case the sample is to be stored below 25°C, the examination has to be started within 6 hours. Exceptionally the sample can be stored at 2 to 8°C (refrigeration) for up to 24 hours.

Store dry at 15°C to 25°C.


If stored under recommended conditions the unopened snap pack has a shelf-life of 3 years after day of production (see expiry date on the label).

Quality Control

Control strains	Inoculum on reference medium	Growth	Colour change to blue-green	Fluorescence	Indole formation
Escherichia coli ATCC® 11775 (WDCM 00090)	10-100	+	+	+	+
Citrobacter freundii ATCC® 8090	10-100	+	+	-	-
Salmonella typhimurium ATCC® 14028 (WDCM 00031)	10-100	+	-	-	-
Klebsiella pneumoniae ATCC® 31488	10-100	+	+	-	-
Pseudomonas aeruginosa ATCC® 10145 (WDCM 00024)	≥1000	No limit	-	-	-

Incubation: 24 h; 35°C; aerobic

Please refer to the actual batch related Certificate of Analysis.

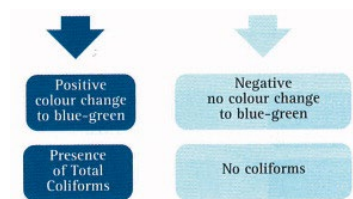

Ordering Information

Product	Cat. No.	Pack size
Readycult® Coliforms 50 (for 50 ml sample)	1.01295 .0001	1 x 20 tests
Readycult® Coliforms 100 (for 100 ml sample)	1.01298 .0001	1 x 20 tests
Bactident® Indole (dropper bottle)	1.11350.0001	1 x 30 ml
CULTURA® Mini-Incubator (100-110 V)	1.15533.0001	1 ea
CULTURA® Mini-Incubator (220-235 V)	1.13311.0001	1 ea
KOVCS Indole Reagent	1.09293.0100	100 ml
UV Lamp (366 nm)	1.13203.0001	1 ea

Merck KGaA

Frankfurter Strasse 250 64293
Darmstadt, Germany Fax: +49
(0) 61 51 / 72-60 80

Find contact information for your
country at:
www.merckmillipore.com/offices

For Technical Service, please visit:
www.merckmillipore.com/techservice

For more information, visit

www.merckmillipore.com/biomonitring

Readycult, Merck, Millipore, and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. Detailed information on trademarks is available via publicly accessible resources.

© 2019 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

The life science business of Merck operates as
MilliporeSigma in the U.S. and Canada.