

QUANTI-Blue™ Solution

Medium for detection and quantification of alkaline phosphatase in standard and HTS assays

Catalog code: rep-qbs, rep-qbs2, rep-qbs3

<https://www.invivogen.com/quant-blue>

For research use only

Version 20C16-MM

PRODUCT INFORMATION

Contents: QUANTI-Blue™ Solution is available in three pack sizes

- **rep-qbs:** 5 x 1 ml of QB reagent and 5 x 1 ml QB buffer, sufficient to prepare QUANTI-Blue™ Solution for **25 x 96-well plates** (500 ml using the standard procedure) or **20 x 1536-well plates** (85 ml using the HTS screening procedure).

- **rep-qbs2:** 10 x 1 ml of QB reagent and 10 x 1 ml QB buffer, sufficient to prepare QUANTI-Blue™ Solution for **50 x 96-well plates** (1 L using the standard procedure) or **40 x 1536-well plates** (170 ml using the HTS screening procedure).

- **rep-qbs3:** 1 x 20 ml bottle of QB reagent and 1 x 20 ml bottle of QB buffer, sufficient to prepare QUANTI-Blue™ Solution for **100 x 96-well plates** (2 L using the standard procedure) or **80 x 1536-well plates** (340 ml using the HTS screening procedure).

Required Material (not provided)

- Sterile water
- Sterile screw cap tube, glass bottle or flask

Storage and stability

- Product is shipped at room temperature. Upon receipt, store QB reagent and QB buffer at -20°C. Product is stable for 1 year at -20°C when properly stored.

- The 20 ml bottles of QB reagent and QB buffer are designed for single use. If required, individual aliquots of QB reagent and QB buffer can be prepared upon receipt or following a single freeze-thaw cycle. Store aliquots at -20°C. **Avoid repeated freeze-thaw cycles.**

Note: During storage, a precipitate may form in the 20 ml bottle of QB reagent. If this occurs, vortex the product until the precipitate disappears. The formation of a precipitate does not affect the activity of the product.

- Reconstituted QUANTI-Blue™ Solution is stable for 2 weeks at 2-8°C and for 2 months at -20°C. Protect QUANTI-Blue™ from light.

Quality Control

Each lot is thoroughly tested to ensure the absence of lot-to-lot variation.

- Physicochemical characterization (including pH, solubility).
- Functional assays using alkaline phosphatase or SEAP-expressing reporter cells.

DESCRIPTION

QUANTI-Blue™ is a colorimetric enzyme assay developed to determine any alkaline phosphatase activity (AP) in a biological sample, such as supernatants of cell cultures. QUANTI-Blue™ Solution changes from pink to a purple-blue color in the presence of AP. Secreted embryonic alkaline phosphatase (SEAP) is a widely used reporter gene. It is a truncated form of placental alkaline phosphatase, a glycosylphosphatidylinositol (GPI)-anchored protein. SEAP is secreted into the cell culture supernatant and therefore offers many advantages over intracellular reporters.

QUANTI-Blue™ is highly sensitive for quantitative measurement. It has a higher saturation threshold than with pNPP (p-nitrophenyl phosphate) resulting in more significant differences between no, low or high AP activity. Another advantage of QUANTI-Blue™ is that it can determine secreted AP activity without disturbing cells, thus allowing the repeated sampling of cell cultures for kinetic studies.

TECHNICAL SUPPORT

InvivoGen USA (Toll-Free): 888-457-5873

InvivoGen USA (International): +1 (858) 457-5873

InvivoGen Europe: +33 (0) 5-62-71-69-39

InvivoGen Asia: +852 3622-3480

E-mail: info@invivogen.com

METHODS

QUANTI-Blue™ Solution has been optimized for use in 96-well plates (standard procedure) and in 1536-well plates (high throughput screening procedure).

A. Standard procedure

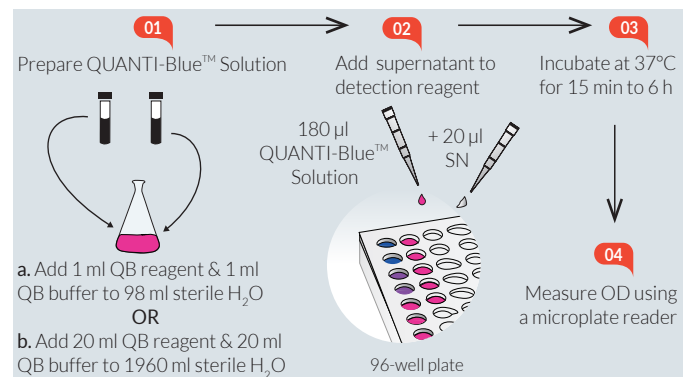


Figure 1. Standard procedure using QUANTI-Blue™ Solution.

The following protocol refers to the use of 96-well plates. Ensure QB reagent and QB buffer are completely thawed before use.

Note: For fast thawing, QB reagent and QB buffer can be placed at 37°C for 2 minutes. Ensure heating at 37°C does **not** exceed 5 minutes.

1. In a sterile bottle or flask, prepare QUANTI-Blue™ Solution by adding:
 - a. 1 ml of QB reagent and 1 ml of QB buffer to 98 ml of sterile water.
 - b. 20 ml of QB reagent and 20 ml of QB buffer to 1960 ml of sterile water.
2. Mix by vortexing and incubate at room temperature for 10 min before use.
3. Use QUANTI-Blue™ Solution immediately or store at 2-8°C or -20°C.
4. Dispense 180 µl of QUANTI-Blue™ Solution per well into a flat-bottom 96-well plate.
5. Add 20 µl of the sample (supernatant of SEAP-expressing cells) or negative control (cell culture medium).
6. Incubate at 37°C for 15 min to 6 h.
7. Measure optical density (OD) at 620-655 nm using a microplate reader.

Note: If the negative control turns purple/blue, it means the fetal bovine serum (FBS) contains alkaline phosphatase. We recommend heating FBS at 56°C for 30 min to inactivate the alkaline phosphatase activity.

For different cell culture plate formats, please refer to the table below:

	96-well plate	24-well plate	12-well plate
QUANTI-Blue™	180 µl	450 µl	900 µl
Supernatant	20 µl	50 µl	100 µl

B. High Throughput Screening (HTS) procedure

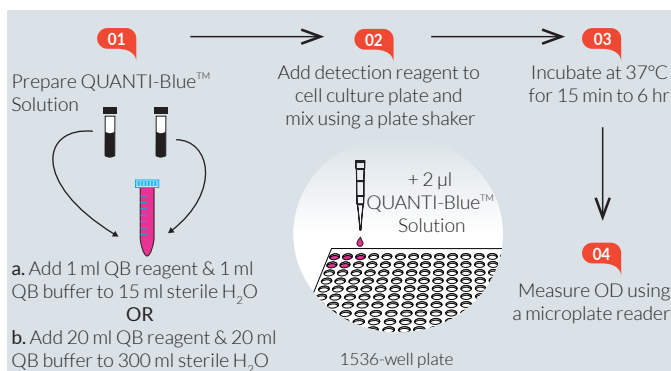


Figure 2. High throughput screening procedure using QUANTI-Blue™ Solution.

This procedure has been optimized for use in HTS screening procedures in 1536-well plates. QUANTI-Blue™ Solution is added directly to the cell suspension to reduce liquid handling.

Ensure QB reagent and QB buffer are completely thawed before use. *Note: For fast thawing, QB reagent and QB buffer can be placed at 37°C for 2 minutes. Ensure heating at 37°C does **not** exceed 5 minutes.*

1. Dispense cell suspension and test compounds into a 1536-well plate in a volume that does not exceed **5 µl** per well. Incubate cells with test compounds for the desired period of time.
2. Prepare QUANTI-Blue™ Solution by adding:
 - a. **1 ml** of QB reagent and **1 ml** of QB buffer to **15 ml** of sterile water in a sterile 50 ml screw cap tube.
 - b. **20 ml** of QB reagent and **20 ml** of QB buffer to **300 ml** of sterile water in a sterile glass bottle or flask.
3. Mix well by vortexing and incubate at room temperature for 10 minutes before use.
4. Use QUANTI-Blue™ Solution immediately or store at 2-8°C or -20°C.
5. Dispense **2 µl** of QUANTI-Blue™ Solution to the wells containing $\leq 5 \mu\text{l}$ of cell culture in a 1536-well plate.
6. Mix using a plate shaker.
7. Incubate at 37°C for 15 min to 6 h.
8. Measure OD at 620-655 nm.

Note: If the negative control turns purple/blue, it means the fetal bovine serum (FBS) contains alkaline phosphatase. We recommend heating FBS at 56°C for 30 min to inactivate the alkaline phosphatase activity.

RELATED PRODUCTS

Product	Catalog Code
pNifTy2-SEAP (Zeo®)	pnifty2-seap
pSELECT-zeo-SEAP	psetz-seap
HEK-Blue™ Detection	hb-det2
Recombinant SEAP Protein	rec-hseap
Reporter cells	
HEK-Blue™ hTLR2	hkb-htlr2
HEK-Blue™ hTLR4	hkb-htlr4
RAW-Blue™ Cells	raw-sp
THP1-Blue™ NF-κB Cells	thp-nfkb
THP1-Blue™ ISG Cells	thp-isg

For a complete list of InvivoGen's Reporter Cell Lines visit <https://www.invivogen.com/reporter-cells>

TECHNICAL SUPPORT

InvivoGen USA (Toll-Free): 888-457-5873
 InvivoGen USA (International): +1 (858) 457-5873
 InvivoGen Europe: +33 (0) 5-62-71-69-39
 InvivoGen Asia: +852 3622-3480
 E-mail: info@invivogen.com